Membranventil

Metall, DN 4 - 50

Diaphragm Valve

Metal, DN 4 - 50

- Ⓓ ORIGINAL EINBAU- UND MONTAGEANLEITUNG
- ⒼB INSTALLATION, OPERATING AND MAINTENANCE INSTRUCTIONS

Membrangröße 8
Diaphragm size 8

Membrangröße 10
Diaphragm size 10

Membrangröße 25 - 50
Diaphragm sizes 25 - 50

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Contents

1	General information	28
2	General safety information	29
2.1	Information for service and operating personnel	29
2.2	Warning notes	29
2.3	Symbols used	30
3	Definition of terms	30
4	Intended area of use	30
5	Technical data	30
6	Order data	32
7	Manufacturer's information	34
7.1	Transport	34
7.2	Delivery and performance	34
7.3	Storage	34
7.4	Tools required	34
8	Functional description	34
9	Construction	34
9.1	Type plate	35
10	Installation and operation	35
10.1	Installing the diaphragm valve	35
10.2	Operation	37
10.3	Setting the seal adjuster	37
10.4	Autoclaving the valve	39
10.4.1	Autoclaving valves with diaphragm sizes 8 - 10	39
10.4.2	Autoclaving valves with diaphragm sizes 25 - 50	39
11	Assembly / disassembly of spare parts	40
11.1	Valve disassembly (removing bonnet from body)	40
11.2	Removing the diaphragm	40
11.3	Mounting the diaphragm	40
11.3.1	General information	40
11.3.2	Mounting a concave diaphragm	41
11.3.3	Mounting a convex diaphragm	42
11.4	Removing the SAK wearing parts kit	42
11.4.1	Removing the 673 8SAK...0P (diaphragm size 8) wearing parts kit	42
11.4.2	Removing the 673 10SAK...1P (diaphragm size 10) wearing parts kit	43
11.4.3	Removing the 673...SAK...P... (diaphragm sizes 25 - 50) wearing parts kit	43

11.5	Mounting the SAK wearing parts kit	44
11.5.1	Mounting the 673 8SAK...0P (diaphragm size 8) wearing parts kit	44
11.5.2	Mounting the 673 10SAK...1P (diaphragm size 10) wearing parts kit	44
11.5.3	Mounting the 673...SAK...P... (diaphragm sizes 25 - 50) wearing parts kit	44
11.6	Bonnet mounting on the valve body	45
12	Commissioning	45
13	Inspection and servicing	45
14	Disassembly	46
15	Disposal	46
16	Returns	46
17	Information	46
18	Troubleshooting / Fault clearance	47
19	Dimensions	48
20	Sectional drawings and spare parts	49
20.1	Diaphragm size 8	49
20.2	Diaphragm size 10	50
20.3	Diaphragm sizes 25 - 50	51
21	EU declaration of conformity	52

1 General information

Prerequisites to ensure that the GEMÜ valve functions correctly:

- x Correct transport and storage
 - x Installation and commissioning by trained personnel
 - x Operation according to these installation, operating and maintenance instructions
 - x Recommended maintenance
- Correct installation, operation, servicing and repair work ensure faultless diaphragm valve operation.

The descriptions and instructions apply to the standard versions. For special versions not described in these installation, operating and maintenance instructions the basic information contained herein applies in combination with any additional special documentation.

All rights including copyright and industrial property rights are expressly reserved.

2 General safety information

The safety information does not take into account:

- x Unexpected incidents and events, which may occur during installation, operation and servicing.
- x Local safety regulations which must be adhered to by the operator and by any additional installation personnel.

2.1 Information for service and operating personnel

The installation, operating and maintenance instructions contain fundamental safety information that must be observed during commissioning, operation and servicing. Non-compliance with these instructions may cause:

- x Personal hazard due to electrical, mechanical and chemical effects.
- x Hazard to nearby equipment.
- x Failure of important functions.
- x Hazard to the environment due to the leakage of dangerous materials.

Prior to commissioning:

- Read the installation, operating and maintenance instructions.
- Provide adequate training for the installation and operating personnel.
- Ensure that the contents of the installation, operating and maintenance instructions have been fully understood by the responsible personnel.
- Define the areas of responsibility.

During operation:

- Keep the installation, operating and maintenance instructions available at the place of use.
- Observe the safety information.
- Use only in accordance with the specifications.
- Any servicing work and repairs not described in the installation, operating

and maintenance instructions must not be performed without consulting the manufacturer first.

⚠ DANGER

Strictly observe the safety data sheets or the safety regulations that are valid for the media used.

In cases of uncertainty:

- x Consult the nearest GEMÜ sales office.

2.2 Warning notes

Wherever possible, warning notes are organised according to the following scheme:

⚠ SIGNAL WORD

Type and source of the danger

- Possible consequences of non-observance.
- Measures for avoiding danger.

Warning notes are always marked with a signal word and sometimes also with a symbol for the specific danger. The following signal words and danger levels are used:

⚠ DANGER

Imminent danger!

- Non-observance will lead to death or severe injury.

⚠ WARNING

Potentially dangerous situation!

- Non-observance can cause death or severe injury.

⚠ CAUTION

Potentially dangerous situation!

- Non-observance can cause moderate to light injury.

CAUTION (WITHOUT SYMBOL)

Potentially dangerous situation!

- Non-observance can cause damage to property.

2.3 Symbols used

	Danger - hot surfaces!
	Danger - corrosive materials!
	Danger - bodily injury!
	Hand: indicates general information and recommendations.
●	Bullet point: indicates the tasks to be performed.
➤	Arrow: indicates the response(s) to tasks.
x	Enumeration sign


3 Definition of terms

Working medium

The medium that flows through the diaphragm valve.

5 Technical data

Working medium	
Corrosive, inert, gaseous and liquid media which have no negative impact on the physical and chemical properties of the body and diaphragm material.	
Temperatures	
Medium temperature	-10 to 100 °C
Sterilisation temperature ⁽¹⁾	
EPDM (code 13/3A)	max. 150 °C ⁽²⁾ , max. 60 min per cycle
EPDM (code 17, 19)	max. 150 °C ⁽²⁾ , max. 180 min per cycle
PTFE/EPDM (code 54)	max. 150 °C ⁽²⁾ , no time limit per cycle
PTFE/EPDM (code 5M)	max. 150 °C ⁽²⁾ , no time limit per cycle
FKM (code 4/4A)	not applicable
¹ The sterilisation temperature is valid for steam (saturated steam) or superheated water.	
² If the sterilisation temperatures listed above are applied to the EPDM diaphragms for longer periods of time, the service life of the diaphragms will be reduced. In these cases, maintenance cycles must be adapted accordingly.	
This also applies to PTFE diaphragms exposed to high temperature fluctuations.	
PTFE diaphragms can also be used as steam barriers; however, this will reduce their service life.	
The maintenance cycles must be adapted accordingly.	
GEMÜ 555 and 505 globe valves are particularly suitable for use in the area of steam generation and distribution.	
The following valve arrangement for interfaces between steam pipes and process pipes has proven itself over time:	
A globe valve for shutting off steam pipes and a diaphragm valve as an interface to the process pipes.	
Ambient temperature	0 to 60 °C

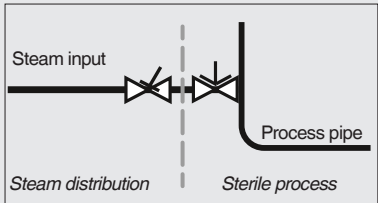4 Intended area of use

- x The GEMÜ 673P9 diaphragm valve is designed for installation in piping systems. It controls a flowing medium by manual operation.
- x **The valve may only be used providing the product technical criteria are complied with (see chapter 5 "Technical data").**
- x Do not paint the bolts and plastic parts of the diaphragm valve!

⚠ WARNING

Use the diaphragm valve only for the intended purpose!

- Otherwise the manufacturer liability and guarantee will be void.
- Use the diaphragm valve only in accordance with the operating conditions specified in the contract documentation and in the installation, operating and maintenance instructions.
- The diaphragm valve must not be used in potentially explosive zones.

Operating pressure [bar]						
		EPDM / FKM		PTFE		
Diaphragm size	Nominal size	Diaphragm material	All valve body materials	Diaphragm material	Forged body	Investment cast or cast iron body
8	DN 4 - 15	3A, 4A, 17, 19	0 - 10	54	0 - 10	0 - 6
10	DN 10 - 20	4, 13, 17, 19	0 - 10	54, 5M	0 - 10	0 - 6
25	DN 15 - 25	4, 13, 17, 19	0 - 10	54, 5M	0 - 10	0 - 6
40	DN 32 - 40	4, 13, 17, 19	0 - 10	54, 5M	0 - 10	0 - 6
50	DN 50 - 65	4, 13, 17, 19	0 - 10	54, 5M	0 - 10	0 - 6

All pressures are gauge pressures. Operating pressure values were determined with static operating pressure applied on one side of a closed valve. Sealing at the valve seat and atmospheric sealing is ensured for the given values. Information on operating pressures applied on both sides and for high purity media on request.

Kv values [m³/h]								
Pipe standard	DIN	EN 10357 series B (formerly DIN 11850 series 1)	EN 10357 series A (formerly DIN 11850 series 2) / DIN 11866 series A	DIN 11850 Series 3	SMS 3008	ASME BPE / DIN 11866 series C	ISO 1127 / EN 10357 series C / DIN 11866 series B	
Connection code	0	16	17	18	37	59	60	
MG	DN							
8	4	0.5	-	-	-	-	-	-
	6	-	-	1.1	-	-	-	1.2
	8	-	-	1.3	-	-	0.6	2.2
	10	-	2.1	2.1	2.1	-	1.3	-
	15	-	-	-	-	-	2.0	-
10	10	-	2.4	2.4	2.4	-	2.2	3.3
	15	3.3	3.8	3.8	3.8	-	2.2	4.0
	20	-	-	-	-	-	3.8	-
25	15	4.1	4.7	4.7	4.7	-	-	7.4
	20	6.3	7.0	7.0	7.0	-	4.4	13.2
	25	13.9	15.0	15.0	15.0	12.6	12.2	16.2
40	32	25.3	27.0	27.0	27.0	26.2	-	30.0
	40	29.3	30.9	30.9	30.9	30.2	29.5	32.8
50	50	46.5	48.4	48.4	48.4	51.7	50.6	55.2
	65	-	-	-	-	62.2	61.8	-

MG = diaphragm size

Kv values determined acc. to DIN EN 60534, inlet pressure 5 bar, Δp 1 bar, stainless steel valve body (forged body) and soft elastomer diaphragm. The Kv values for other product configurations (e.g. other diaphragm or body materials) may differ. In general, all diaphragms are subject to the influences of pressure, temperature, the process and their tightening torques. Therefore the Kv values may exceed the tolerance limits of the standard.

The Kv value curve (Kv value dependent on valve stroke) can vary depending on the diaphragm material and duration of use.

6 Order data

Body configuration	Code
Tank bottom valve body	B**
2/2-way body	D
T body	T*
* For dimensions see T Valves brochure	
** Dimensions and versions on request	

Connection	Code
Butt weld spigots	
Spigots DIN	0
Spigots EN 10357 series B (formerly DIN 11850 series 1)	16
Spigot EN 10357 series A (formerly DIN 11850 series 2) / DIN 11866 series A	17
Spigots DIN 11850 series 3	18
Spigots JIS-G 3447	35
Spigots JIS-G 3459	36
Spigots SMS 3008	37
Spigots BS 4825 Part 1	55
Spigot ASME BPE / DIN 11866 series C	59
Spigot ISO 1127 / EN 10357 series C / DIN 11866 series B	60
Spigots ANSI/ASME B36.19M Schedule 10s	63
Spigots ANSI/ASME B36.19M Schedule 5s	64
Spigots ANSI/ASME B36.19M Schedule 40s	65
Threaded connections	
Threaded sockets DIN ISO 228	1
Threaded spigots DIN 11851	6
Cone spigot and union nut DIN 11851	6K
Aseptic unions on request	
Flanges	
Flanges EN 1092 / PN16 / form B, length EN 558, series 1, ISO 5752, basic series 1	8
Flanges ANSI Class 150 RF, length MSS SP-88	38
Flanges ANSI Class 125/150 RF, length EN 558, series 1, ISO 5752, basic series 1	39
Clamp connections	
Clamps ASME BPE for pipe ASME BPE, length ASME BPE	80
Clamps DIN 32676 series B for pipe EN ISO 1127, length EN 558, series 7	82
Clamps ASME BPE for pipe ASME BPE, length EN 558, series 7	88
Clamps DIN 32676 series A for pipe DIN 11850, length EN 558, series 7	8A
Clamps SMS 3017 for pipe SMS 3008, length EN 558, series 7	8E
Clamps DIN 32676 series C, length FTF ASME BPE	8P
Clamps DIN 32676 series C, length FTF EN 558 series 7	8T
Aseptic clamps on request	

Valve body material	Code
EN-GJS-400-18-LT (SG iron 40.3) PFA lined	17
EN-GJS-400-18-LT (SG iron 40.3) PP lined	18
1.4435, investment casting	C3
1.4408, investment casting	37
1.4435 (316L), forged body	40
1.4435 (BN2), forged body $\Delta Fe < 0.5\%$	42
1.4539, forged body	F4

Diaphragm material	Code
FKM	4 4A*
EPDM	13 3A*
EPDM	17 17*
EPDM	19
EPDM	36
PTFE/EPDM, one-piece	54*
PTFE/EPDM, two-piece	5M
* for diaphragm size 8	
Material complies with FDA requirements, except codes 4 and 4A	

Control function	Code
Manually operated	0

Bonnet version	Code
With seal adjuster, black handwheel diaphragm size 8	0P9
With seal adjuster, black handwheel diaphragm size 10	1P9
With seal adjuster, black handwheel diaphragm size 25	2P9
With seal adjuster, black handwheel diaphragm size 40	3P9
With seal adjuster, black handwheel diaphragm size 50	4P9

Internal surface finishes for forged and block material bodies ¹

Readings for Process Contact Surfaces	Mechanically polished ²		Electropolished	
	Hygienic class DIN 11866	Code	Hygienic class DIN 11866	Code
Ra ≤ 0.80 µm	H3	1502	HE3	1503
Ra ≤ 0.60 µm	-	1507	-	1508
Ra ≤ 0.40 µm	H4	1536	HE4	1537
Ra ≤ 0.25 µm ³	H5	1527	HE5	1516

Readings for Process Contact Surfaces acc. to ASME BPE 2016 ⁴	Mechanically polished ²		Electropolished	
	ASME BPE Surface Designation	Code	ASME BPE Surface Designation	Code
Ra Max. = 0.76 µm (30 µinch)	SF3	SF3	-	-
Ra Max. = 0.64 µm (25 µinch)	SF2	SF2	SF6	SF6
Ra Max. = 0.51 µm (20 µinch)	SF1	SF1	SF5	SF5
Ra Max. = 0.38 µm (15 µinch)	-	-	SF4	SF4

Internal surface finishes for investment cast bodies

Readings for Process Contact Surfaces	Mechanically polished ²	
	Hygienic class DIN 11866	Code
Ra ≤ 6.30 µm	-	1500
Ra ≤ 0.80 µm	H3	1502
Ra ≤ 0.60 µm ⁵	-	1507

¹ Surface finishes of customized valve bodies may be limited in special cases.

² Or any other finishing method that meets the Ra value (acc. to ASME BPE).

³ The smallest possible Ra finish for pipe connections with an internal pipe diameter < 6 mm is 0.38 µm.

⁴ When using these surfaces, the bodies are marked according to the specifications of ASME BPE.

The surfaces are only available for valve bodies which are made of materials (e.g. GEMÜ material codes 40, 41, F4, 44) and use connections (e.g. GEMÜ connection codes 59, 80, 88) according to ASME BPE.

⁵ Not possible for GEMÜ connection code 59, DN 8 and GEMÜ connection code 0, DN 4.

Ra acc. to DIN EN ISO 4288 and ASME B46.1

Order example	673	8	D	60	40	17	0	0P9	1537
Type	673								
Nominal size		8							
Body configuration (code)			D						
Connection (code)				60					
Valve body material (code)					40				
Diaphragm material (code)						17			
Control function (code)							0		
Bonnet version (code)								0P9	
Surface finish (code)									1537

7 Manufacturer's information

7.1 Transport

- Only transport the diaphragm valve by suitable means. Do not drop. Handle carefully.
- Dispose of packing material according to relevant local or national disposal regulations / environmental protection laws.

7.2 Delivery and performance

- Check that all parts are present and check for any damage immediately upon receipt.
- The scope of delivery is apparent from the dispatch documents and the design from the order number.
- The performance of the diaphragm valve is checked at the factory.

7.3 Storage

- Store the diaphragm valve free from dust and moisture in its original packaging.
- Store the diaphragm valve in "open" position.
- Avoid UV rays and direct sunlight.
- Maximum storage temperature: 40 °C.
- Solvents, chemicals, acids, fuels or similar fluids must not be stored in the same room as valves and their spare parts.

7.4 Tools required

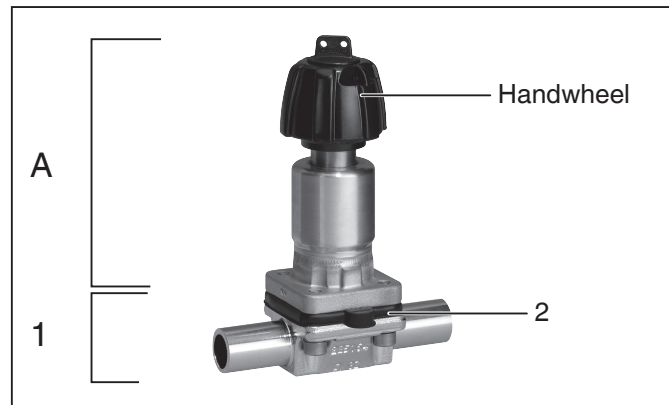
- The tools required for installation and assembly are **not** included in the scope of delivery.
- Use appropriate, functional and safe tools.

8 Functional description

GEMÜ 673P9 is a metal diaphragm valve with a 2/2-way, T or tank bottom valve body or in multi-port design. The manually operated GEMÜ 673P9 diaphragm valve has better tightness characteristics due to


its unique seal system. It has a temperature resistant plastic handwheel. Bonnet housing and internals are made all of stainless steel. An integral optical position indicator and a seal adjuster to increase service life of the diaphragm are included as standard. The valve body and the diaphragm are available in various designs as shown in the data sheet. The valve has CIP / SIP / COP, sterilizing and autoclave capabilities.

9 Construction


Construction - Diaphragm size 8

1	Valve body
2	Diaphragm
A	Bonnet

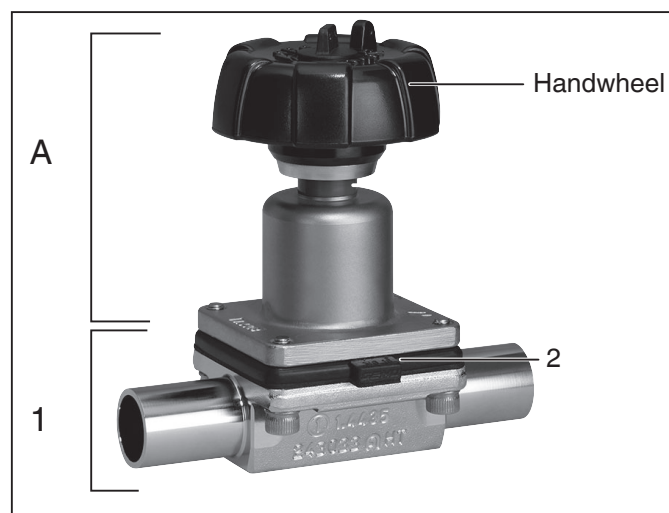Construction - Diaphragm size 10

1	Valve body
2	Diaphragm
A	Bonnet

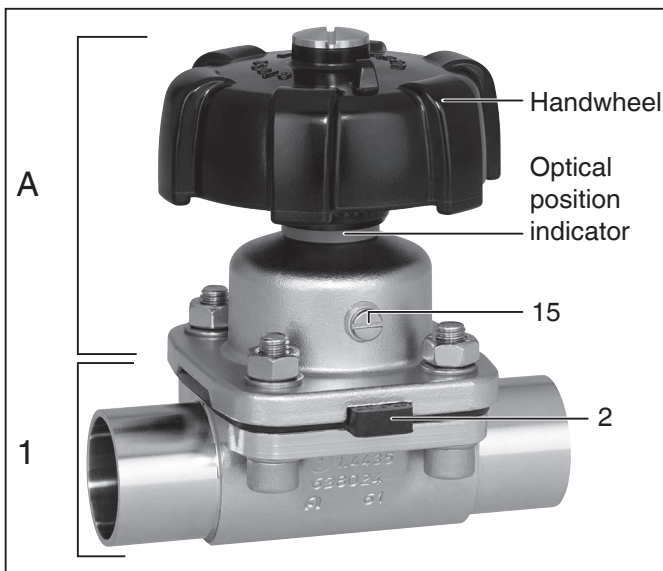Construction - Diaphragm sizes 25-50

1	Valve body
2	Diaphragm
15	Special screw
A	Bonnet

9.1 Type plate

Device version	Design in accordance with order data		Device-specific data	
	673	8D60405400P9	PS 10,0 bar	Year of manufacture
	1502			
	ERC DE	2020		
	88701322	12103529	10001	
Item number	Traceability number		Serial number	

The month of manufacture is encoded in the traceability number and can be obtained from GEMÜ.

The product was manufactured in Germany.

10 Installation and operation

Prior to installation:

- Ensure that valve body and diaphragm material are appropriate and compatible to handle the working medium.
- **Check the suitability prior to the installation.**
See chapter 5 "Technical data".

10.1 Installing the diaphragm valve

⚠ WARNING

The equipment is subject to pressure!

- Risk of severe injury or death!
- Only work on depressurized plant.

⚠ WARNING

Corrosive chemicals!

- Risk of caustic burns!
- Wear appropriate protective gear when installing.

⚠ CAUTION

Hot plant components!

- Risk of burns!
- Only work on plant that has cooled down.

⚠ CAUTION

Never use the valve as a step or an aid for climbing!

- This entails the risk of slipping-off or damaging the valve.

CAUTION

Do not exceed the maximum permissible pressure!

- Take precautionary measures to avoid possible pressure surges (water hammer).

- Installation work must only be performed by trained personnel.
- Use appropriate protective gear as specified in plant operator's guidelines.

Installation location:

⚠ CAUTION

- Do not apply external force to the valve.
- Choose the installation location so that the valve cannot be used as a foothold (climbing aid).
- Lay the pipeline so that the valve body is protected against transverse and bending forces, and also vibrations and tension.
- Only install the valve between matching aligned pipes.

- x Direction of the working medium: optional.
- x Mounting position of the diaphragm valve: optional.

Installation:

1. Ensure the suitability of the valve for each respective use. The valve must be appropriate for the piping system operating conditions (medium, medium concentration, temperature and pressure) and the prevailing ambient conditions. Check the technical data of the valve and the materials.
2. Shut off plant or plant component.
3. Secure against recommissioning.
4. Depressurize the plant or plant component.
5. Completely drain the plant (or plant component) and let it cool down until the temperature is below the media vaporization temperature and scalding can be ruled out.
6. Correctly decontaminate, rinse and ventilate the plant or plant component.

Installation - Butt weld spigots:

1. Adhere to good welding practices!
2. Disassemble the bonnet with the diaphragm before welding the valve body into the pipeline (see chapter 11.1).
3. Allow butt weld spigots to cool down.
4. Reassemble the valve body and the bonnet with diaphragm (see chapter 11.6).

Installation - Clamp connections:

- When installing clamp connections, insert a gasket between the body clamp and the adjacent piping clamp and join them using the appropriate clamp fitting. The gasket and the clamp for clamp connections are not included in the scope of delivery.

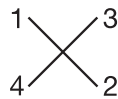
Important:

Butt weld spigots / Clamp connections:

Angle of rotation for welding into pipeline to enable optimised draining see brochure "Angle of rotation for 2/2-way valve bodies" (on request or www.gemu-group.com).

Installation - Flanges:

1. Pay attention to clean, undamaged sealing surfaces on the mating flanges.
2. Align flanges carefully before installing them.
3. Centre the seals accurately.
4. Connect the valve flange and the piping flange using appropriate sealing material and matching bolting. Sealing material and bolts are not included in the scope of delivery.
5. Use all flange holes.
6. Only use connector elements made of approved materials!
7. Tighten the bolts diagonally!

Installation - Threaded connections:

- Screw the threaded connections into the piping in accordance with valid standards.
- Screw the diaphragm valve body into the piping, use appropriate thread sealant. The thread sealant is not included in the scope of delivery.

Observe appropriate regulations for connections!

After the installation:

- Reactivate all safety and protective devices.

10.2 Operation

Optical position indicator:

Diaphragm size 8

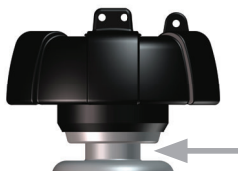
Valve open

Valve closed

Diaphragm size 10

Valve open

Valve closed

Diaphragm sizes 25 - 50

Valve open

Valve closed

⚠ CAUTION

Handwheel can become hot during operation!

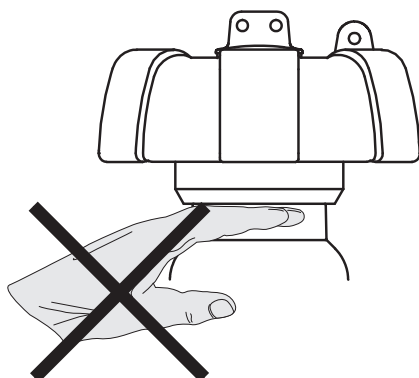
- Risk of burns!
- Ensure protective gloves are worn when operating handwheel.

⚠ CAUTION

Rising handwheel!

- Danger of crushing fingers.

10.3 Setting the seal adjuster

Important:

Only set the seal adjuster when the valve is completely assembled (with diaphragm and valve body) and in a cold condition!

The valves type GEMÜ 673P9 have a mechanical seal adjuster as standard to protect the sealing diaphragm.

Standard setting:

The valve is sealed when the handwheel is completely closed (turned fully clockwise).

Setting the seal adjuster:

Procedure:

1. Open the valve approx. 50 %.
2. Unscrew special screw 6.

Diaphragm size 8

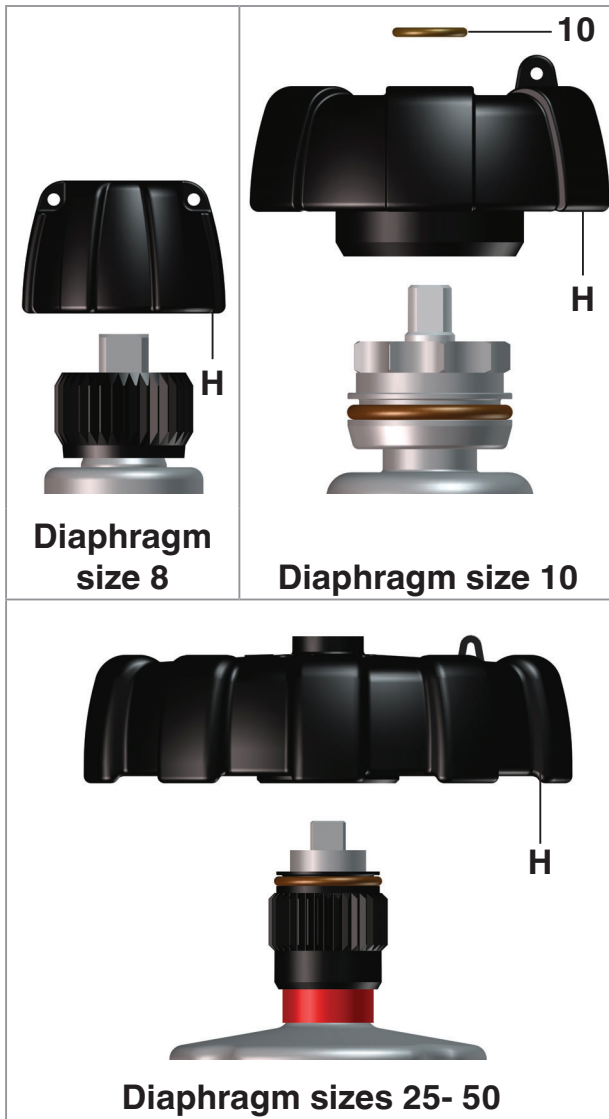
Diaphragm size 10


Diaphragm sizes 25- 50


3. Pull up handwheel **H** and O-ring **10** if existing.

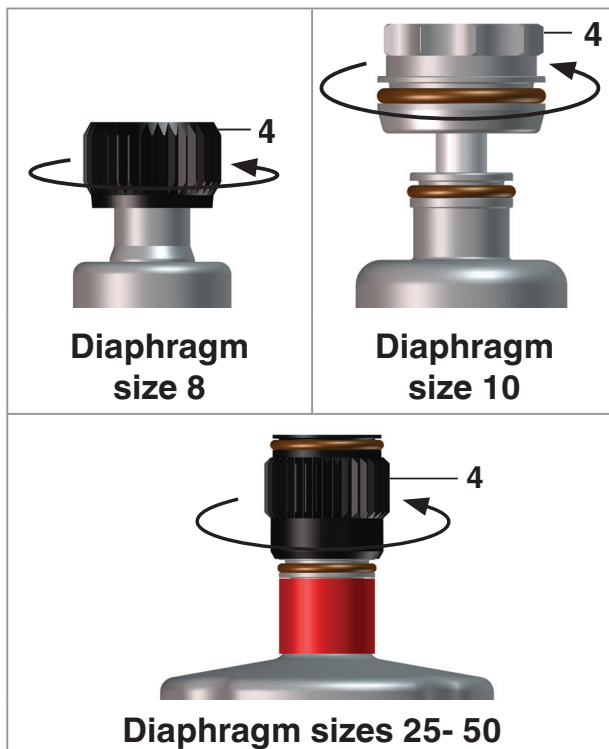
4. Unscrew distance piece **4**.

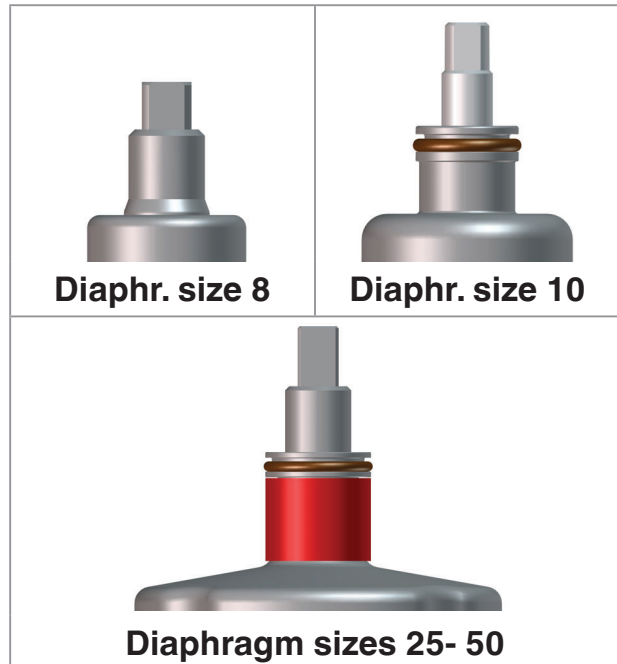5. Dependent on the operating conditions, regrease the threaded spindle, especially a valve that is autoclaved. GEMÜ recommends the grease Boss-Fluorine Y 108/00 (99099484).

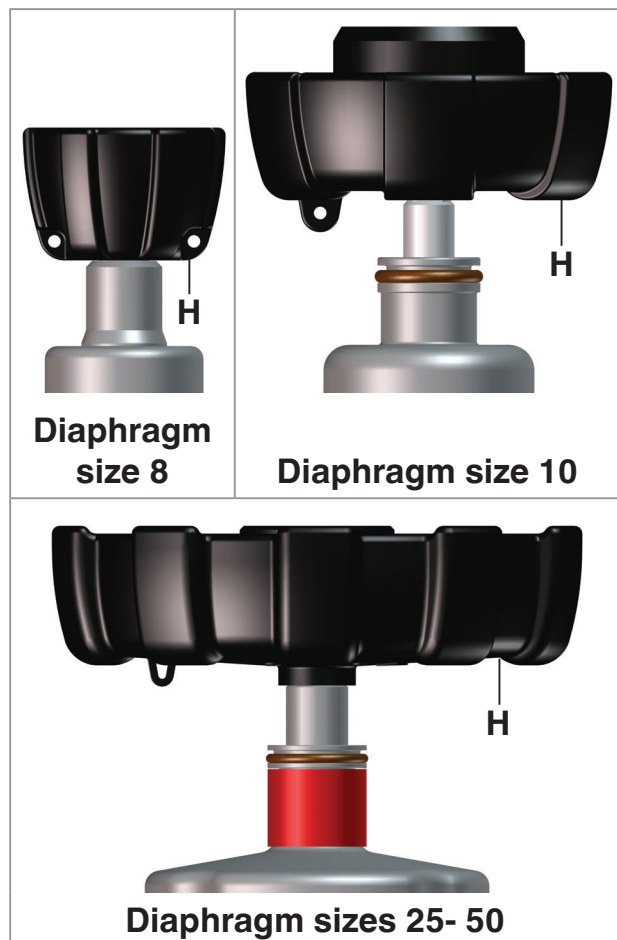
6. Place the handwheel **H** upside down on the square of the threaded spindle.


7. Close the valve gently by turning handwheel **H** ("CLOSED").

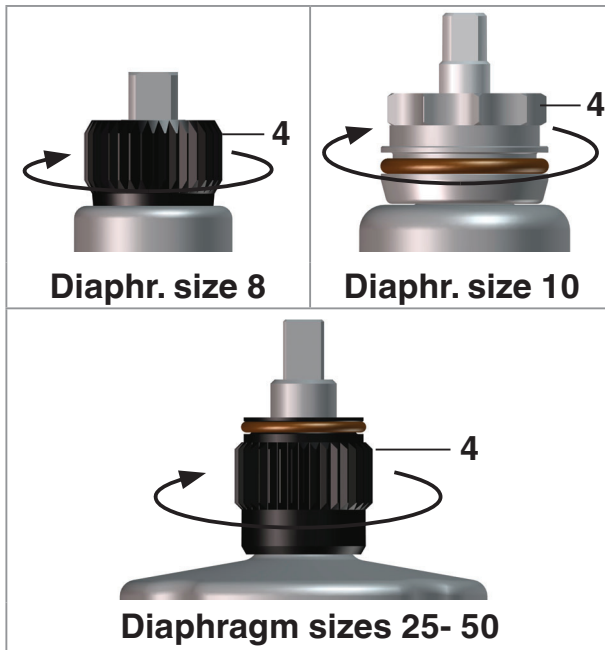8. Pull off handwheel **H** from the threaded spindle.
9. Position distance piece **4** (bevel downwards) on the threaded spindle and screw it down clockwise until it sits on the valve bonnet.

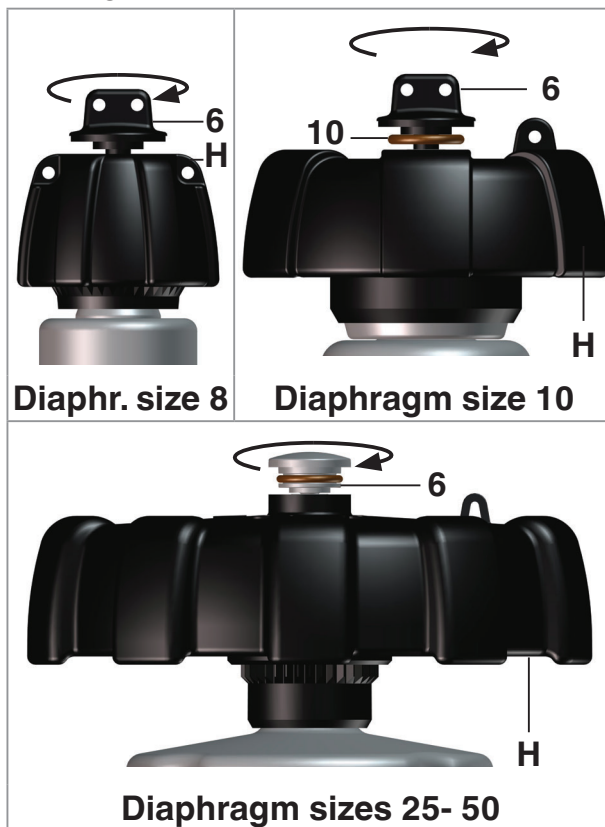10. Push handwheel **H** in its original position on the square of the threaded spindle (observe the teeth of the adjusting ring **4** and the handwheel **H**). Tighten it with special screw **6** (insert O-ring **10** if existing).

10.4 Autoclaving the valve

⚠ CAUTION

High temperatures during autoclaving!

➤ Damage to seals.

- Move the bonnet to the open position.
- Before autoclaving ensure that special screw **15** is tightened hand tight (only for diaphragm sizes 25 - 50).

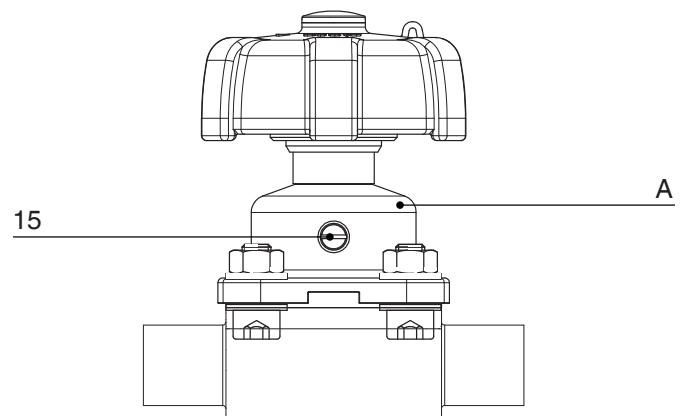
Diaphragm sizes 8 and 10 have no leak detection hole.

10.4.1 Autoclaving valves with diaphragm sizes 8 - 10

1. Move bonnet **A** to the open position.
2. Autoclave valve **A**.

10.4.2 Autoclaving valves with diaphragm sizes 25 - 50


1. Move bonnet **A** to the open position.
2. Hand tighten special screw **15** without tools.
3. Autoclave valve **A**.
4. Loosen special screw **15** after autoclaving.

Only loosen special screw **15** where this is necessary for the further work process.

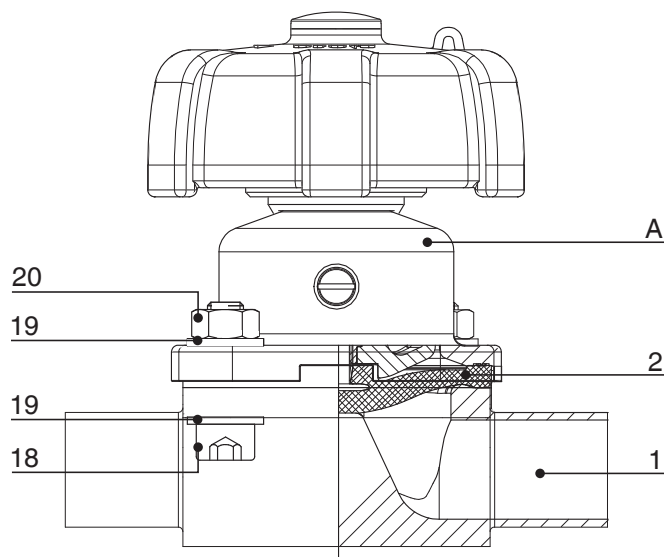11 Assembly / disassembly of spare parts

11.1 Valve disassembly (removing bonnet from body)

1. Move bonnet **A** to the open position.
2. Remove bonnet **A** from valve body **1**.
3. Move bonnet **A** to the closed position.

Important:

After disassembly, clean all parts of contamination (do not damage parts). Check parts for potential damage, replace if necessary (only use genuine parts from GEMÜ).

11.2 Removing the diaphragm

Important:

Before removing the diaphragm, please remove the bonnet, see "Valve disassembly (removing bonnet from body)".

1. Unscrew the diaphragm or pull it out (diaphragm size 8).
2. Clean all parts of the remains of product and contamination. Do not scratch or damage parts during cleaning!
3. Check all parts for potential damage.
4. Replace damaged parts (only use genuine parts from GEMÜ).

11.3 Mounting the diaphragm

11.3.1 General information

Important:

Mount the correct diaphragm that suits the valve (suitable for medium, medium concentration, temperature and pressure). The diaphragm is a wearing part. Check the technical condition and function of the diaphragm valve before commissioning and during the whole term of use. Carry out checks regularly and determine the check intervals in accordance with the conditions of use and / or the regulatory codes and provisions applicable for this application.

Important

(diaphragm sizes 10 - 50):

If the diaphragm is not screwed into the adapter far enough, the closing force is transmitted directly onto the diaphragm pin and not via the compressor. This will cause damage and early failure of the diaphragm and thus leakage of the valve. If the diaphragm is screwed in too far no perfect sealing at the valve seat will be achieved. The function of the valve is no longer ensured.

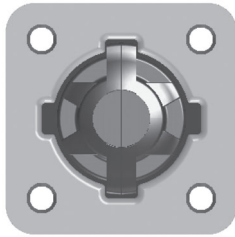
Important:

Incorrectly mounted diaphragm may cause valve leakage / emission of medium. In this case remove the diaphragm, check the complete valve and diaphragm and reassemble again proceeding as described above.

The compressor is fixed to the spindle for all diaphragm sizes.


Diaphragm size 8:

Compressor and bonnet flange seen from below:


Diaphragm size 10:

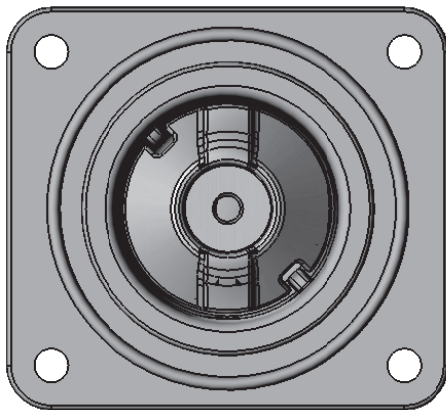Compressor and bonnet flange seen from below:

Diaphragm sizes 25 - 50:

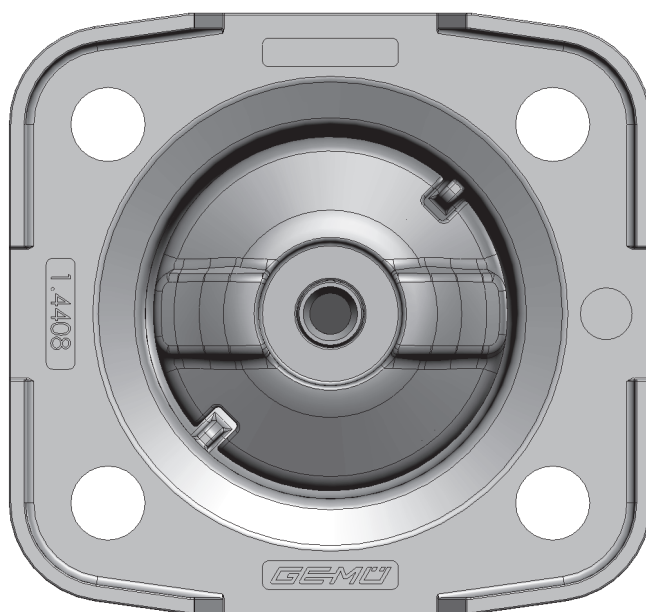
Compressor and bonnet flange seen from below:

11.3.2 Mounting a concave diaphragm

Diaphragm size 8

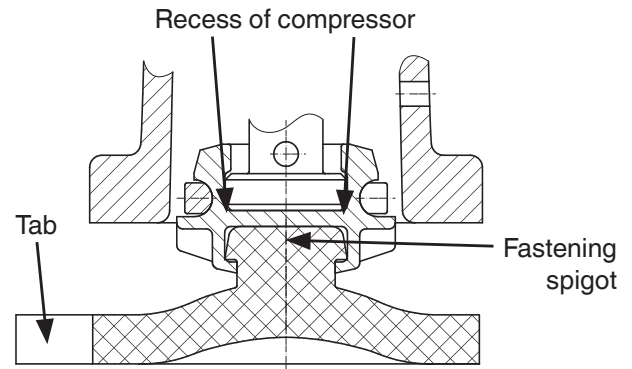Push-fit diaphragm:

1. Move bonnet **A** to the closed position.
2. Place the diaphragm **2** with the fastening spigot in an inclined position at the recess of the compressor.

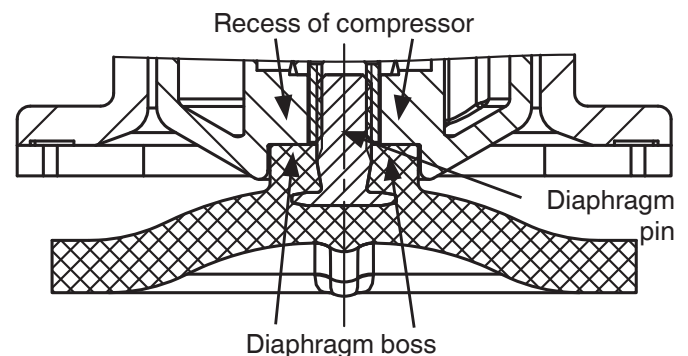
Important:

Do not use any greases or lubricants!

3. Turn the diaphragm as manual force is applied to push the spigot into the compressor.
4. Align diaphragm tab (identifying manufacturer and material) in parallel to compressor weir.

Diaphragm sizes 10 - 50

Threaded pin type diaphragm:


1. Move bonnet **A** to the closed position.
2. Screw new diaphragm tightly into the compressor manually.
3. Check if the diaphragm boss fits closely in the recess of the compressor.
4. If it is difficult to screw it in, check the thread, replace damaged parts (only use genuine parts from GEMÜ).
5. When clear resistance is felt turn back the diaphragm anticlockwise until its bolt


holes are in correct alignment with the bolt holes of the bonnet.

11.3.3 Mounting a convex diaphragm

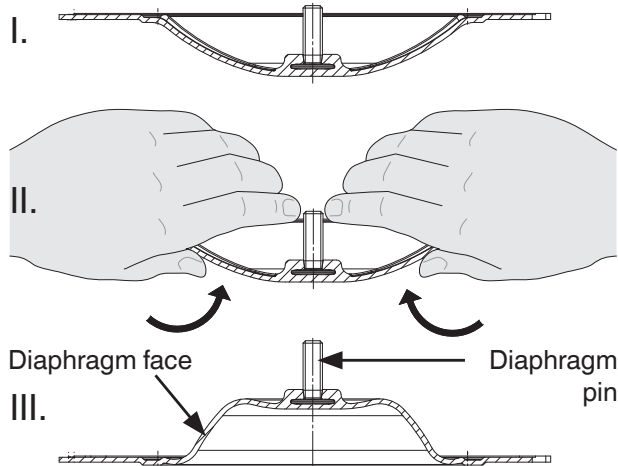1. Move bonnet **A** to the closed position.
2. Invert the new diaphragm face manually; use a clean, padded mat with bigger nominal sizes.

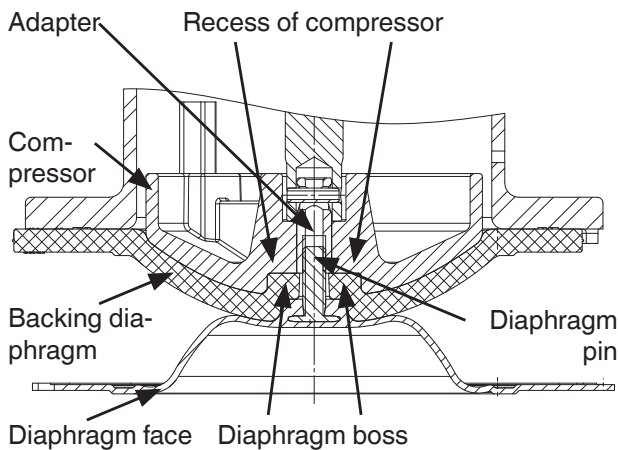3. Position the new backing diaphragm onto the compressor.
4. Position the diaphragm face onto the backing diaphragm.
5. Screw diaphragm face tightly into the compressor manually. The diaphragm boss must fit closely in the recess of the compressor.

6. If it is difficult to screw it in, check the thread, replace damaged parts.
7. When clear resistance is felt turn back the diaphragm anticlockwise until its bolt holes are in correct alignment with the bolt holes of the bonnet.
8. Press the diaphragm face tightly onto the backing diaphragm manually so that it returns to its original shape and fits closely on the backing diaphragm.

11.4 Removing the SAK wearing parts kit

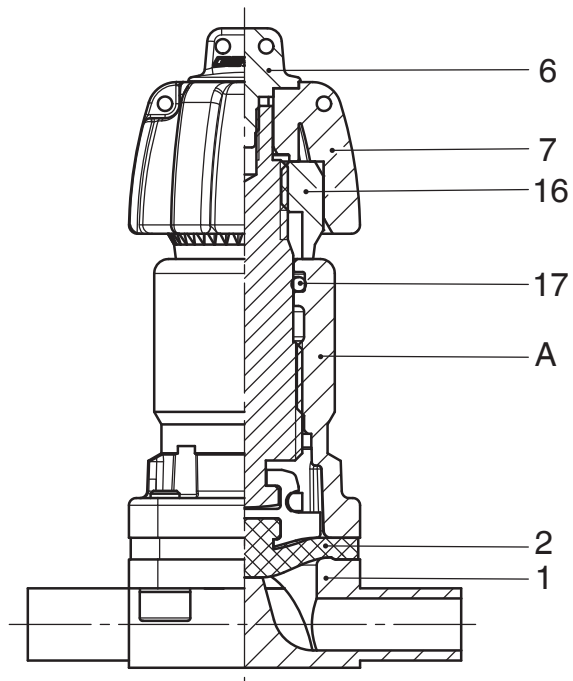
For information on the accessories for the appropriate wearing parts kit, see chapter 20 "Sectional drawings and spare parts".

11.4.1 Removing the 673 8SAK...0P (diaphragm size 8) wearing parts kit


Before removing the wearing parts kit, remove the bonnet, see "Valve disassembly (removing bonnet from body)".

1. Unscrew the special screw **6**.
2. Pull the handwheel **7** upwards from the threaded spindle.
3. Unscrew the distance piece **16** from the threaded spindle.
4. Remove the bonnet **A** from the valve body **1** (see chapter 11.1 "Valve disassembly" (removing bonnet from body)).
5. Pull out and remove the diaphragm **2** (see chapter 11.2 "Removing the diaphragm").
6. Place the handwheel upside down on the square of the threaded spindle.
7. Turn the handwheel **7** clockwise.
 - The threaded spindle is unscrewed from the bonnet housing **A**.


8. Remove the O-ring **17** from the bonnet **A** using a suitable tool.
9. Clean all parts of the remains of product and contamination. Do not scratch or damage parts during cleaning!
10. Check all parts for potential damage.
11. Replace damaged parts (only use genuine parts from GEMÜ).

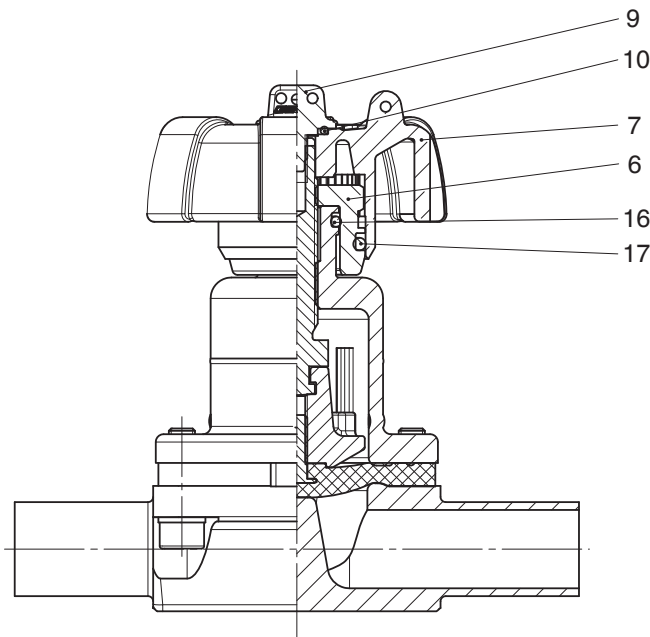11.4.2 Removing the 673 10SAK...1P (diaphragm size 10) wearing parts kit

1. Unscrew the special screw **9**.
2. Remove the O-ring **10** from the handwheel **7**.
3. Pull the handwheel **7** upwards from the threaded spindle.
4. Unscrew the distance piece **6** from the threaded spindle.
5. Remove the O-ring **17** from the distance piece **6**.
6. Remove the O-ring **16** from the bonnet.
7. Clean all parts of the remains of product and contamination. Do not scratch or damage parts during cleaning!
8. Check all parts for potential damage.
9. Replace damaged parts (only use genuine parts from GEMÜ).

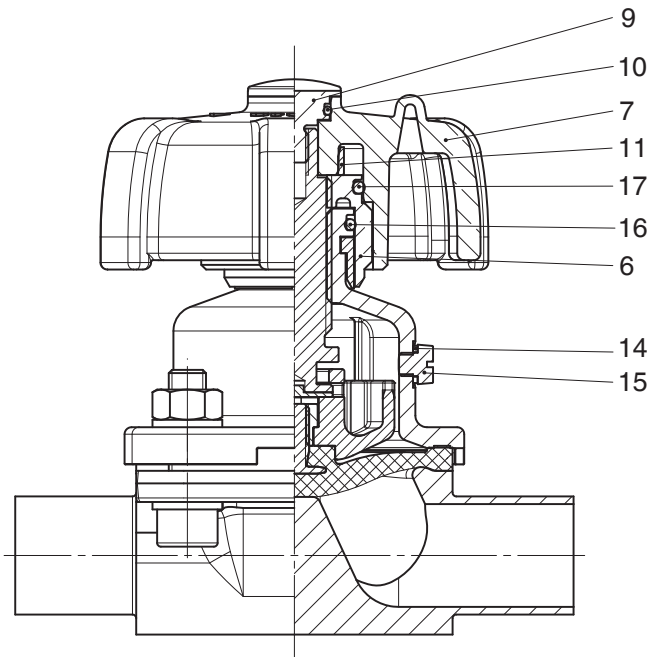11.4.3 Removing the 673...SAK...P... (diaphragm sizes 25 - 50) wearing parts kit

1. Unscrew the special screw **9**.
2. Remove the O-ring **10** from the special screw **9**.
3. Pull the handwheel **7** upwards from the threaded spindle.

The clamping ring **11** is already pre-fitted in the handwheel **7**.

4. Unscrew the distance piece **6** from the threaded spindle.
5. Remove the O-ring **17** from the distance piece **6**.
6. Remove the O-ring **16** from the bonnet.
7. Unscrew the special screw **15** using a flathead screwdriver.
8. Remove the gasket **14**.
9. Clean all parts of the remains of product and contamination. Do not scratch or damage parts during cleaning!
10. Check all parts for potential damage.
11. Replace damaged parts (only use genuine parts from GEMÜ).

11.5 Mounting the SAK wearing parts kit

11.5.1 Mounting the 673 8SAK...0P (diaphragm size 8) wearing parts kit

	See sectional drawing in chapter 11.4.1 "Removing the 673 8SAK...0P (diaphragm size 8) wearing parts kit".
--	--

1. Remove the wearing parts kit (see chapter 11.4.1 "Removing the 673 8SAK...0P (diaphragm size 8) wearing parts kit").
2. Move bonnet **A** to the closed position.
3. Fit the O-ring **17** in the bonnet using a suitable tool.
4. Screw / push the new diaphragm tightly into the compressor manually (see chapter 11.3 "Mounting the diaphragm").
5. Connect the bonnet **A** to the valve body **1** (see chapter 11.6 "Mounting the bonnet on the valve body").
6. Turn the threaded spindle into the bonnet housing.
7. Screw the distance piece **16** onto the threaded spindle.
8. Place the handwheel **7** on the square of the threaded spindle.
9. Fasten the handwheel **7** using the special screw **6**.
10. If necessary, reset the seal adjuster (see chapter 10.3 "Setting the seal adjuster").

11.5.2 Mounting the 673 10SAK...1P (diaphragm size 10) wearing parts kit

	See sectional drawing in chapter 11.4.2 "Removing the 673 10SAK...1P (diaphragm size 10) wearing parts kit".
--	--

1. Remove the wearing parts kit (see chapter 11.4.2 "Removing the 673 10SAK...1P (diaphragm size 10) wearing parts kit").
2. Move bonnet **A** to the closed position.
3. Place the O-ring **16** on the bonnet.
4. Fit the O-ring **17** on the distance piece **6**.

5. Screw the distance piece **6** onto the threaded spindle.
6. Place the handwheel **7** on the square of the threaded spindle.
7. Place the O-ring **10** on the handwheel **7**.
8. Fasten the handwheel **7** using the special screw **9**.
9. If necessary, reset the seal adjuster (see chapter 10.3 "Setting the seal adjuster").

11.5.3 Mounting the 673...SAK...P... (diaphragm sizes 25 - 50) wearing parts kit

	See sectional drawing in chapter 11.4.3 "Removing the 673...SAK...P... (diaphragm size 25-50) wearing parts kit".
---	---


1. Remove the wearing parts kit (see chapter 11.4.3 "Removing the 673...SAK...P... (diaphragm size 25-50) wearing parts kit").
2. Move bonnet **A** to the closed position.
3. Place the O-ring **16** on the bonnet.
4. Fit the O-ring **17** on the distance piece **6**.
5. Screw the distance piece **6** onto the threaded spindle.
6. Place the clamping ring **11** in the handwheel **7**.
7. Place the handwheel **7** on the square of the threaded spindle.

	The clamping ring 11 is already pre-fitted in the handwheel 7 .
---	---

8. Fit the O-ring **10** on the special screw **9**.
9. Fasten the handwheel **7** using the special screw **9**.
10. Insert the gasket **14**.
11. Tighten the special screw **15** using a flathead screwdriver.
12. If necessary, reset the seal adjuster (see chapter 10.3 "Setting the seal adjuster").

11.6 Bonnet mounting on the valve body

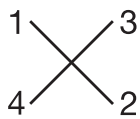1. Move bonnet **A** to the closed position.
2. Open bonnet **A** approx. 50 %.
3. Clean all parts of the remains of product and contamination. Do not scratch or damage parts during cleaning!
4. Position bonnet **A** with the mounted diaphragm **2** on the valve body **1**, take care to align the diaphragm weir and valve body weir.
5. Tighten bolts **18**, washers **19** and nuts **20** by hand (hand tight only).
6. Fully tighten the bolts **18** with nuts **20** diagonally.

7. Ensure that the diaphragm **2** is compressed evenly (approx. 10 - 15 %, visible by an even bulge to the outside).
8. Check tightness of completely assembled valve.

Important:

Service and maintenance:
Diaphragms degrade in the course of time. After valve disassembly / assembly check that the bolts **18** and nuts **20** on the body are tight and retighten as necessary (at the very latest after the first sterilisation process).

⚠ CAUTION

Protect against leakage!

- Provide precautionary measures against exceeding the maximum permitted pressures caused by pressure surges (water hammer).

Prior to cleaning or commissioning the plant:

- Check the tightness and the function of the diaphragm valve (close and reopen the diaphragm valve).
- If the plant is new and after repairs rinse the piping system with a fully opened diaphragm valve (to remove any harmful foreign matter).

Cleaning:

- x The plant operator is responsible for selecting the cleaning material and performing the procedure.

Important:

Service and maintenance:
Diaphragms degrade in the course of time. After valve disassembly / assembly check that the bolts **18** and nuts **20** on the body are tight and retighten as necessary (at the very latest after the first sterilisation process).

12 Commissioning

⚠ WARNING

Corrosive chemicals!

- Risk of caustic burns!
- Check the tightness of the media connections prior to commissioning!
- Use only the appropriate protective gear when performing the tightness check.

13 Inspection and servicing

⚠ WARNING

The equipment is subject to pressure!

- Risk of severe injury or death!
- Only work on depressurized plant.

⚠ CAUTION

Hot plant components!

- Risk of burns!
- Only work on plant that has cooled down.

⚠ CAUTION

- Servicing and maintenance work must only be performed by trained personnel.
- GEMÜ shall assume no liability whatsoever for damages caused by improper handling or third-party actions.
- In case of doubt, contact GEMÜ before commissioning.

1. Use appropriate protective gear as specified in plant operator's guidelines.
2. Shut off plant or plant component.
3. Secure against recommissioning.
4. Depressurize the plant or plant component.

The operator must carry out regular visual examination of the valves dependent on the operating conditions and the potential danger in order to prevent leakage and damage. The valve also has to be disassembled in corresponding intervals and checked for wear (see chapter 11 "Assembly / disassembly of spare parts").

Important:

Service and maintenance:
Dependent on the operating conditions, regrease the threaded spindle, especially a valve that is autoclaved.
See chapter 10.3, point 5.

14 Disassembly

Disassembly is performed observing the same precautionary measures as for installation.

- Disassemble the diaphragm valve (see chapter 11.1 "Valve disassembly (removing bonnet from body)").

15 Disposal

- All valve parts must be disposed of according to relevant local or national disposal regulations / environmental protection laws.
- Pay attention to adhered residual material and gas diffusion from penetrated media.

16 Returns

- Clean the diaphragm valve.
- Request a goods return declaration form from GEMÜ.
- Returns must be made with a completed goods return declaration.

If not completed, GEMÜ cannot process

x credits or

x repair work

but will dispose of the goods at the operator's expense.

Note for returns:

Legal regulations for the protection of the environment and personnel require that the completed and signed goods return declaration is included with the dispatch documents. Returned goods can be processed only when this declaration is completed.

17 Information

Note on staff training:

Please contact us at the address on the last page for staff training information.

Should there be any doubts or misunderstandings in the preceding text, the German version of this document is the authoritative document!


18 Troubleshooting / Fault clearance

Fault	Possible cause	Fault clearance
Medium escapes from leak detection hole* (leak detection hole only existing for diaphragm size 25 - 50, if the special screw 15 was opened or removed)	Shut off diaphragm faulty	Check shut off diaphragm for damage, replace diaphragm if necessary
Valve doesn't open or doesn't open fully	Bonnet faulty	Replace bonnet
	Shut off diaphragm incorrectly mounted	Remove bonnet, check diaphragm mounting, replace if necessary
Valve leaks downstream (doesn't close or doesn't close fully)	Operating pressure too high	Operate valve with operating pressure specified in data sheet
	Foreign matter between shut off diaphragm and valve body weir	Remove bonnet, remove foreign matter, check shut off diaphragm and valve body weir for damage and replace if necessary
	Valve body weir leaking or damaged	Check valve body weir for damage, if necessary replace valve body
	Shut off diaphragm faulty	Check shut off diaphragm for damage, replace diaphragm if necessary
	Seal adjuster is incorrectly set	Readjust the seal adjuster
Valve leaks between bonnet and valve body	Shut off diaphragm incorrectly mounted	Remove bonnet, check diaphragm mounting, replace if necessary
	Bolting between valve body and bonnet loose	Retighten bolting between valve body and bonnet
	Shut off diaphragm faulty	Check shut off diaphragm for damage, replace diaphragm if necessary
	Valve body / bonnet damaged	Replace valve body / bonnet
Valve body connection to piping leaks	Incorrect installation	Check installation of valve body in piping
	Bolting / threaded connections loose	Tighten bolting / threaded connections
	Sealing material faulty	Replace sealing material
Valve body leaks	Valve body faulty or corroded	Check valve body for damage, replace valve body if necessary
Handwheel cannot be turned	Bonnet faulty	Replace bonnet
	Threaded spindle seized	Dependent on the operating conditions, regrease the threaded spindle, especially a valve that is autoclaved; replace bonnet if necessary. See chapter 10.3, point 5.


* see chapter 20 "Sectional drawings and spare parts"

19 Dimensions

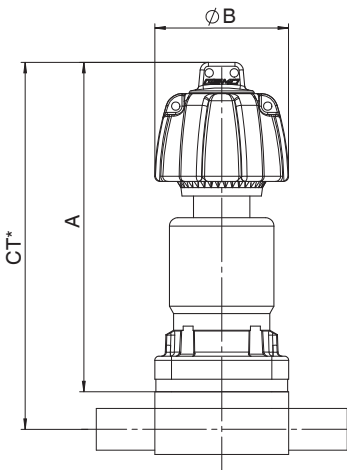
Diaphragm size 8

Diaphragm size 10

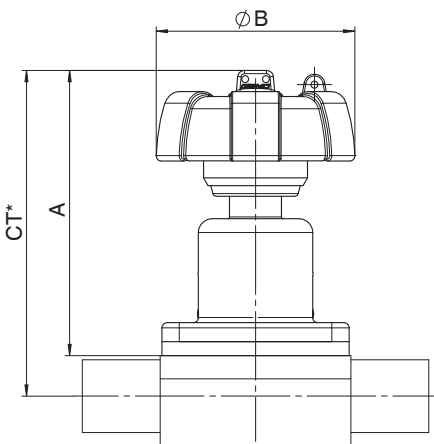Diaphragm size 25 - 50

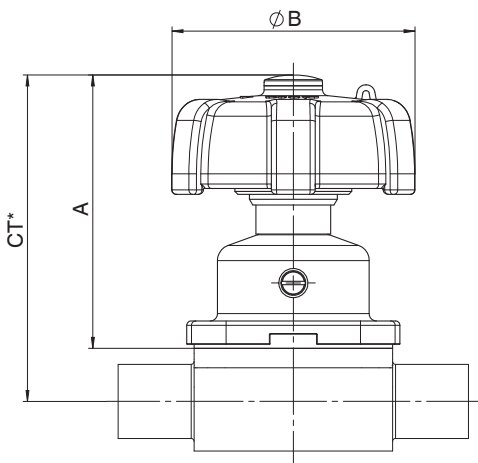
* CT = A + H1 (see body dimensions in data sheet)

Bonnet dimensions			
Diaphragm size	A	Ø B	Weight [kg]
8	80	32	0.20
10	83	60	0.30
25	94	90	0.60
40	121	114	1.25
50	131	140	2.00


20 Sectional drawings and spare parts

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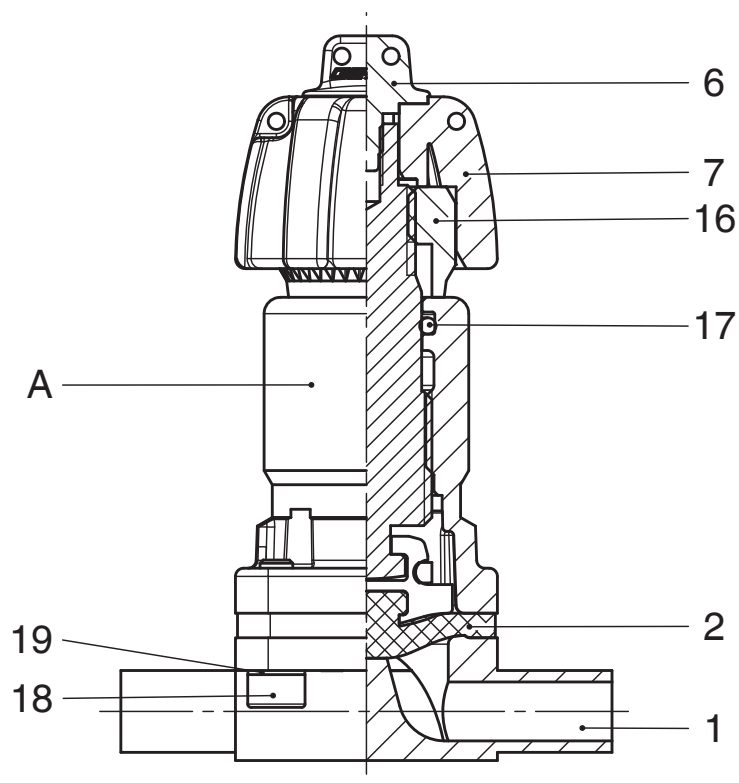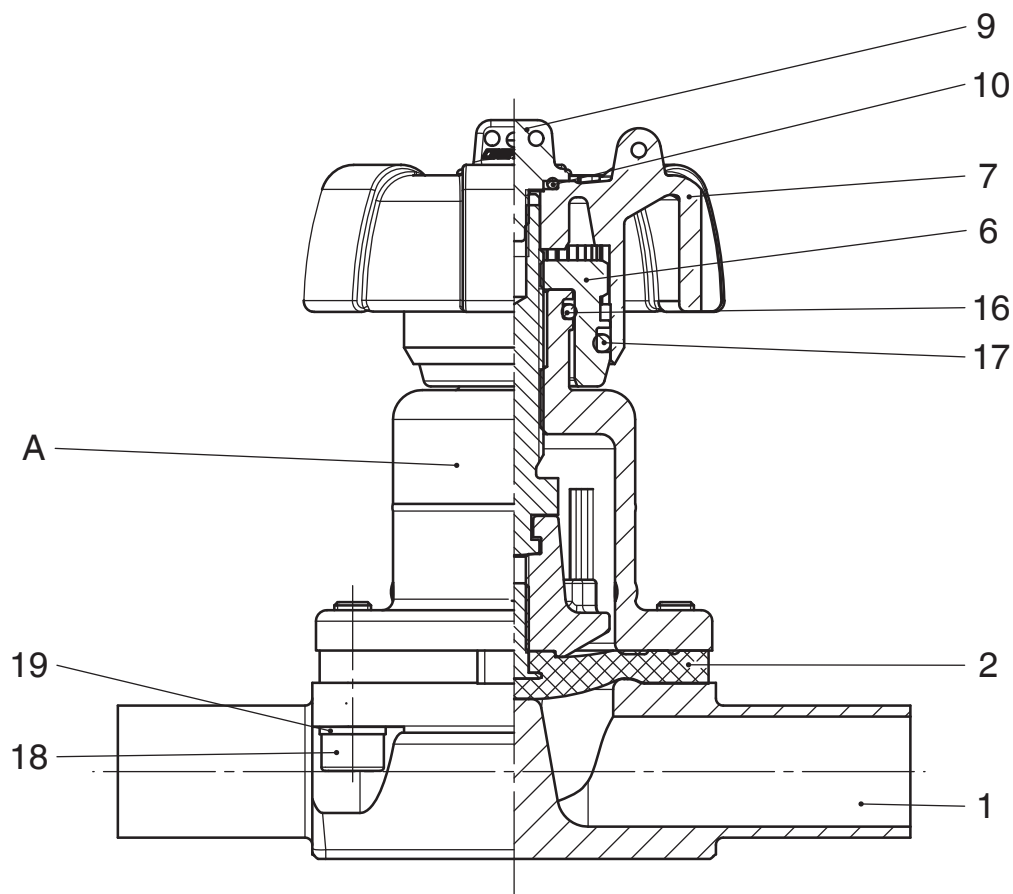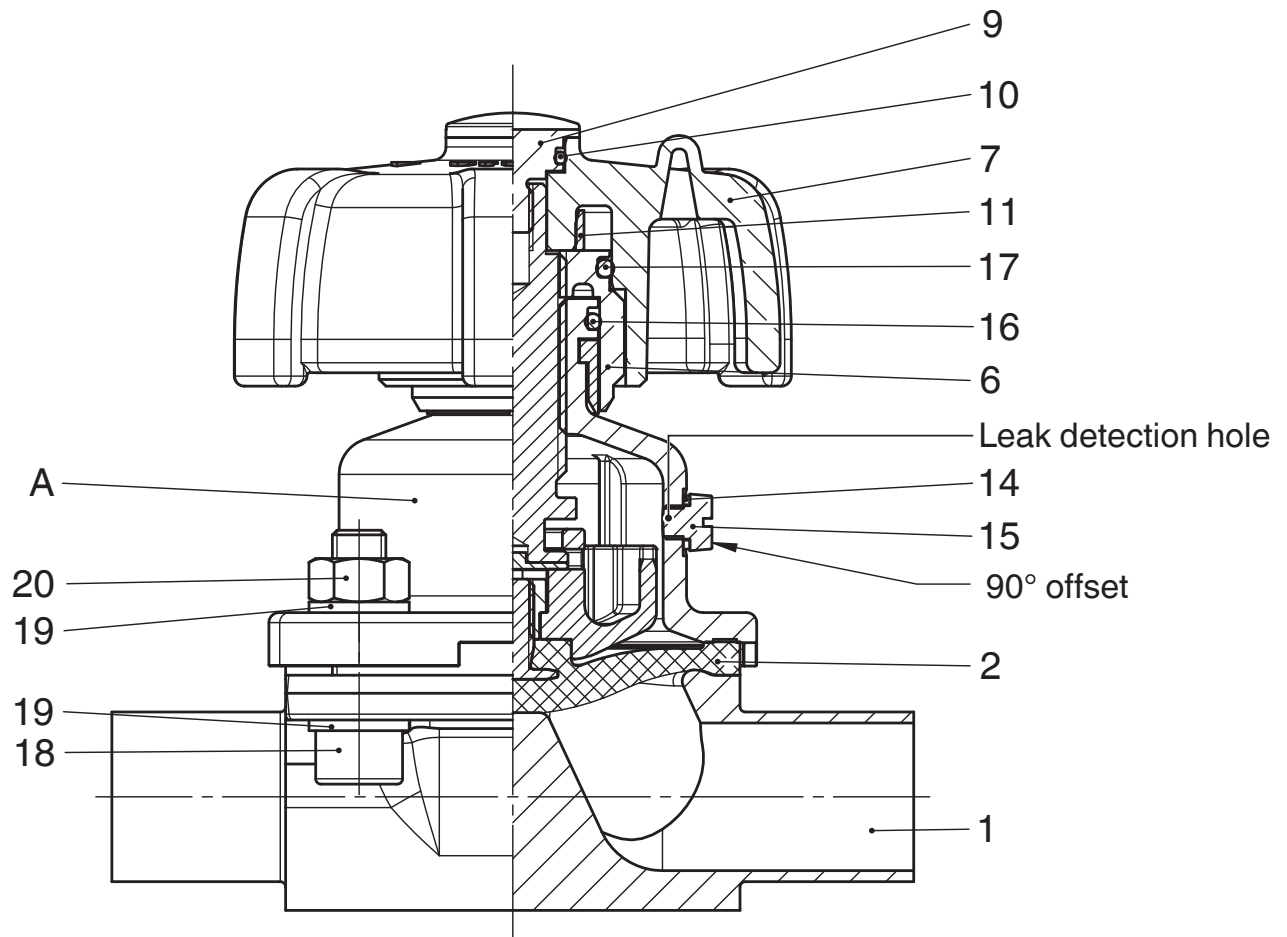
declare that the equipment listed below complies with the safety requirements of the Pressure Equipment Directive 2014/68/EU.

Description of the equipment - product type

Diaphragm valve
GEMÜ 673P9

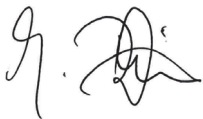
Notified body: TÜV Rheinland Industrie Service GmbH
Number: 0035
Certificate no.: 01 202 926/Q-02 0036
Applied standards: AD 2000

Conformity assessment procedure:
Module H1

Note for equipment with a nominal size \leq DN 25:

The products are developed and produced according to GEMÜ process instructions and quality standards which comply with the requirements of ISO 9001 and of ISO 14001.

According to section 4, paragraph 3 of the Pressure Equipment Directive 2014/68/EU these products must not be identified by a CE-label.

Joachim Brien
Head of Technical Department

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Hereby we, **GEMÜ Gebr. Müller Apparatebau GmbH & Co. KG**