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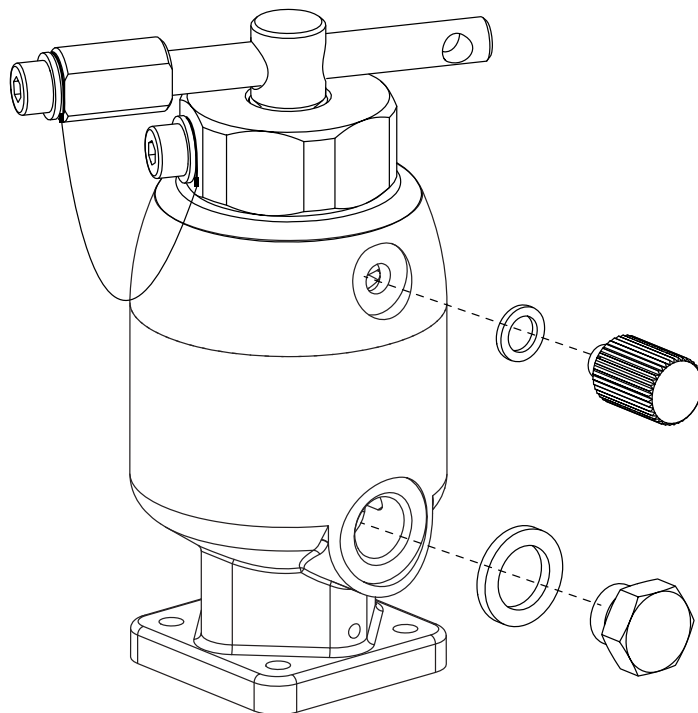
**Special version GEMÜ 650
with an M5 adapter as an exhaust air duct possibility
Metal, diaphragm sizes 8**

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GB ADDITION TO INSTALLATION, OPERATING AND
MAINTENANCE INSTRUCTIONS

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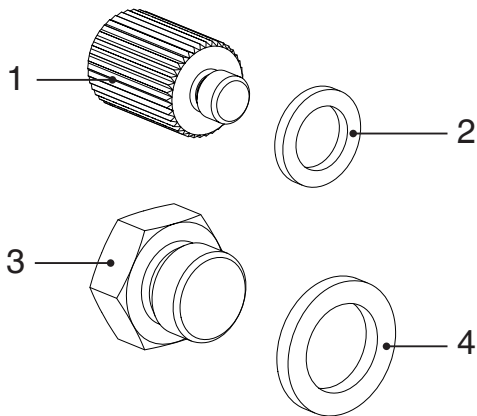
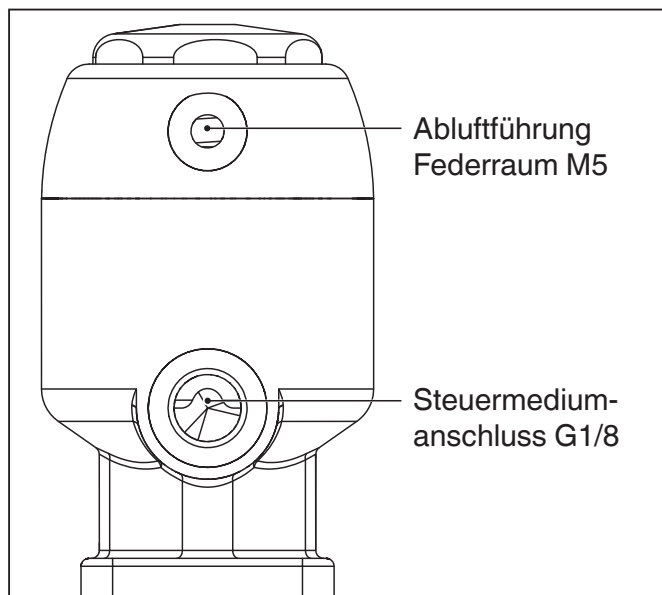
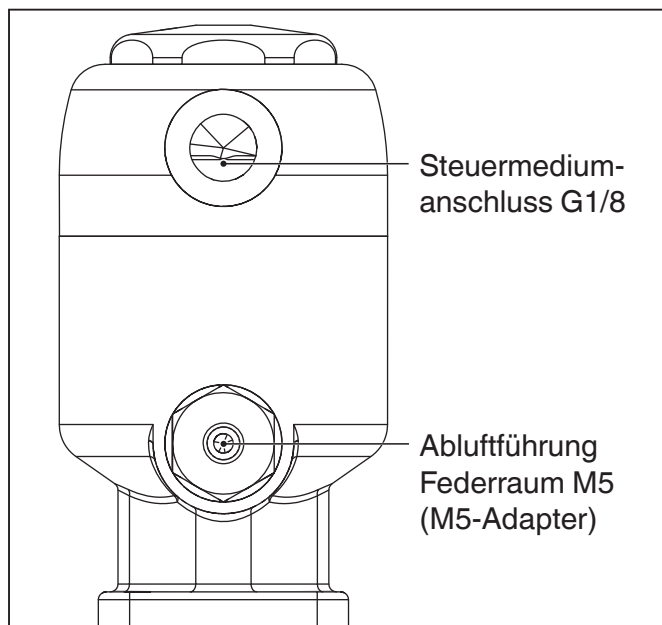
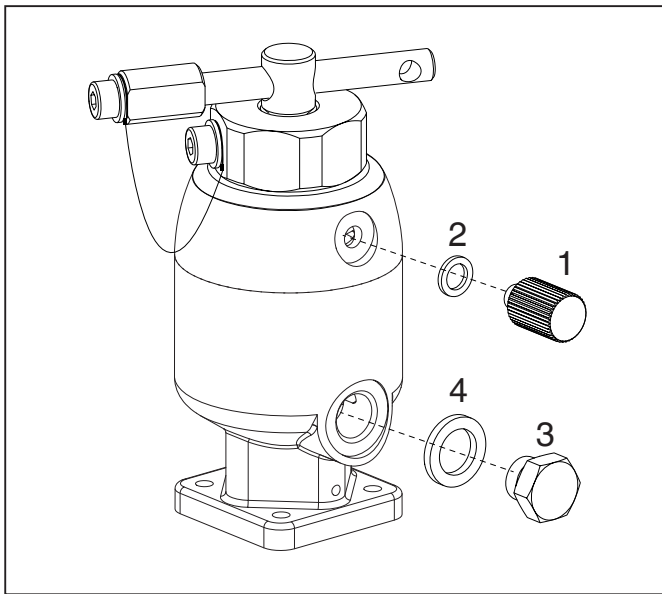
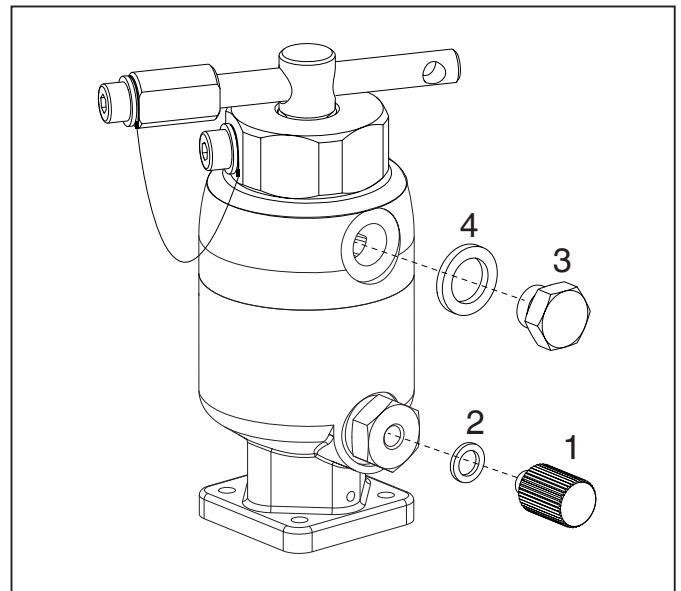
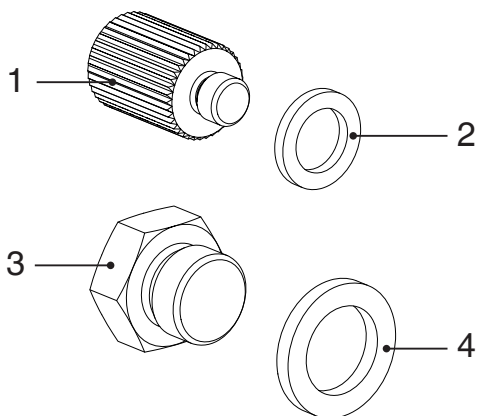
1 General information

Observe the installation, operating and maintenance instructions of the GEMÜ 650.

In addition to their normal function, diaphragm valves with K-no. 5759 allow the exhaust air produced in the spring chamber and in the chamber behind the diaphragm to be ducted via an M5 adapter.

2 Autoclaving kit


The following is included in the scope of delivery:

- x 1x Special screw 1
- x 1x Gasket 2
- x 1x Threaded plug 3
- x 1x Gasket 4


3 Connectors

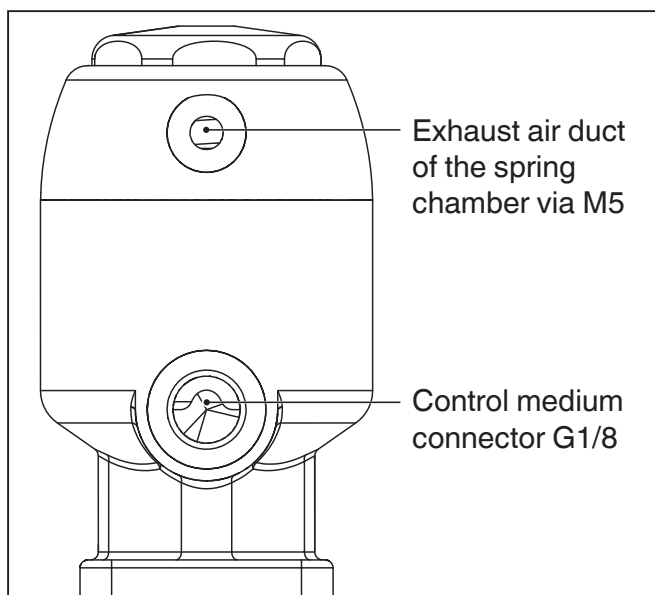3.1 Control function 1

GEMÜ 650, MG 8, control function 1

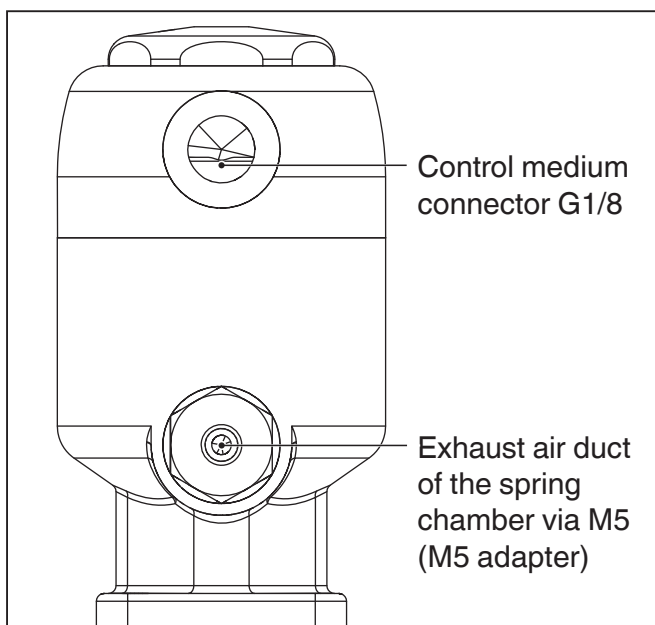
3.2 Control function 2

GEMÜ 650, MG 8, control function 2

4 Autoclaving the valve

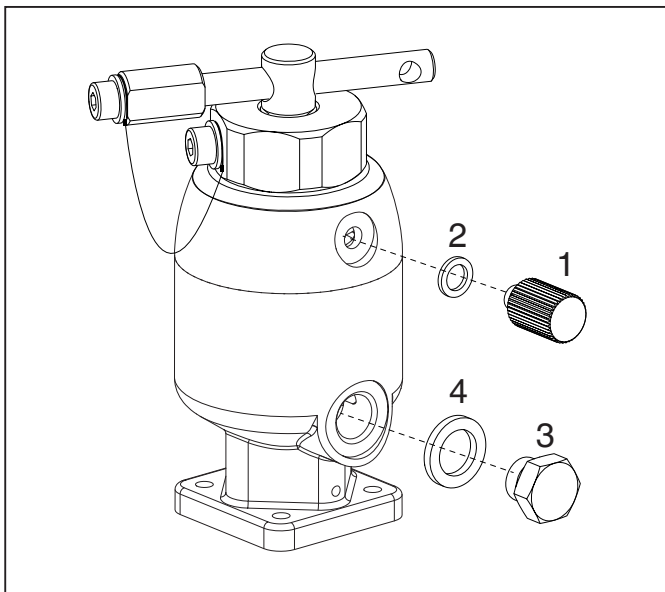
4.1 Control function 1

GEMÜ 650, MG 8, control function 1

1. Move actuator to the open position.
2. Mount the tool to keep actuator open GEMÜ 1107 (see separate assembly instructions).
3. Mount gasket **2** and special screw **1** to exhaust air duct.
4. Mount gasket **4** and threaded plug **3** to control medium connector.
5. Autoclave the valve.
6. Remove gasket **2** and special screw **1** from exhaust air duct.
7. Remove gasket **4** and threaded plug **3** from control medium connector.

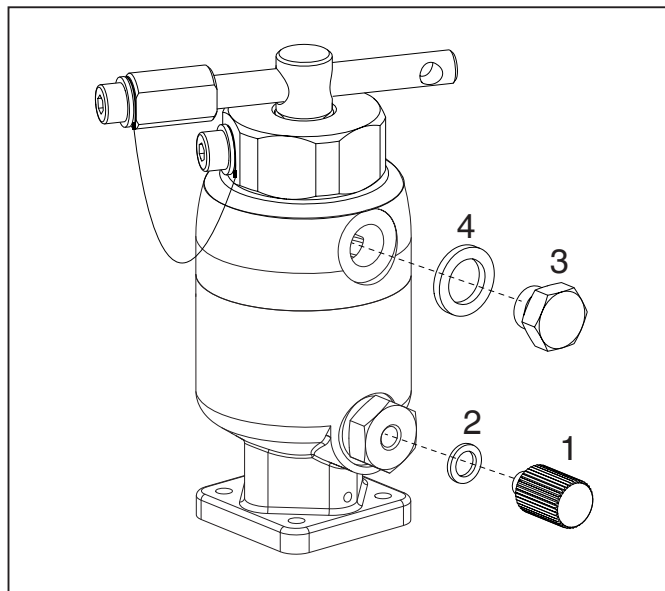4.2 Control function 2

GEMÜ 650, MG 8, control function 2

1. Mount the tool to keep actuator open GEMÜ 1107 (see separate assembly instructions).
2. Mount gasket **2** and special screw **1** to exhaust air duct.
3. Mount gasket **4** and threaded plug **3** to control medium connector.
4. Autoclave the valve.
5. Remove gasket **2** and special screw **1** from exhaust air duct.
6. Remove gasket **4** and threaded plug **3** from control medium connector.