## Membranventil

Metall, DN 4 - 100

## Diaphragm valve

Metal, DN 4 - 100

- Ⓓ ORIGINAL EINBAU- UND MONTAGEANLEITUNG
- Ⓖ INSTALLATION, OPERATING AND MAINTENANCE INSTRUCTIONS

DN 100 "T"

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Actuator version "T"

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Actuator version "D"

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## Contents

<b>1</b>	<b>General information</b>	<b>27</b>
<b>2</b>	<b>General safety information</b>	<b>27</b>
2.1	Information for service and operating personnel	28
2.2	Warning notes	28
2.3	Symbols used	28
<b>3</b>	<b>Definition of terms</b>	<b>29</b>
<b>4</b>	<b>Intended area of use</b>	<b>29</b>
<b>5</b>	<b>Technical data</b>	<b>29</b>
<b>6</b>	<b>Order data (2/2-way valves)</b>	<b>33</b>
<b>7</b>	<b>Manufacturer's information</b>	<b>35</b>
7.1	Transport	35
7.2	Delivery and performance	35
7.3	Storage	35
7.4	Tools required	35
<b>8</b>	<b>Functional description</b>	<b>35</b>
<b>9</b>	<b>Construction</b>	<b>35</b>
9.1	Product label	36
<b>10</b>	<b>Installation and operation</b>	<b>36</b>
10.1	Installing the diaphragm valve	36
10.2	Control functions	37
10.3	Connecting the control medium	38
10.4	Optical position indicator	38
<b>11</b>	<b>Fitting/removing spare parts</b>	<b>39</b>
11.1	Valve disassembly (removing actuator from body)	39
11.2	Removing the diaphragm	39
11.3	Mounting the diaphragm	39
11.3.1	General information	39
11.3.2	Mounting a concave diaphragm	41
11.3.3	Mounting a convex diaphragm	42
11.4	Actuator mounting on the valve body	42
<b>12</b>	<b>Commissioning</b>	<b>43</b>
<b>13</b>	<b>Inspection and maintenance</b>	<b>43</b>
<b>14</b>	<b>Cleaning and sterilization</b>	<b>44</b>
<b>15</b>	<b>Disassembly</b>	<b>44</b>
<b>16</b>	<b>Disposal</b>	<b>44</b>
<b>17</b>	<b>Returns</b>	<b>44</b>
<b>18</b>	<b>Information</b>	<b>44</b>
<b>19</b>	<b>Troubleshooting / Fault clearance</b>	<b>45</b>
<b>20</b>	<b>Sectional drawings and spare parts</b>	<b>46</b>
<b>21</b>	<b>Declaration of Incorporation</b>	<b>50</b>
<b>22</b>	<b>EU Declaration of Conformity</b>	<b>51</b>

## 1 General information

Prerequisites to ensure that the GEMÜ valve functions correctly:

- x Correct transport and storage
- x Installation and commissioning by trained personnel
- x Operation according to these installation, operating and maintenance instructions
- x Recommended maintenance

Correct installation, operation, maintenance and repair work ensure faultless diaphragm valve operation.

The descriptions and instructions apply to the standard versions. For special versions not described in these installation, operating and maintenance instructions the basic information contained herein applies in combination with any additional special documentation.

All rights including copyrights or industrial property rights are expressly reserved.

## 2 General safety information

The safety information does not take into account:

- x Unexpected incidents and events, which may occur during installation, operation and maintenance.
- x Local safety regulations which must be adhered to by the operator and by any additional installation personnel.

2.1 Information for service and operating personnel

The installation, operating and maintenance instructions contain fundamental safety information that must be observed during commissioning, operation and maintenance. Non-compliance with these instructions may cause:

- x Personal hazard due to electrical, mechanical and chemical effects.
- x Hazard to nearby equipment.
- x Failure of important functions.
- x Hazard to the environment due to the leakage of dangerous materials.

Prior to commissioning:

- Read the installation, operating and maintenance instructions.
- Provide adequate training for installation and operating personnel.
- Ensure that the contents of the installation, operating and maintenance instructions have been fully understood by the responsible personnel.
- Define the areas of responsibility.

During operation:

- Keep the installation, operating and maintenance instructions available at the place of use.
- Observe the safety information.
- Use only in accordance with the specifications.
- Any maintenance work and repairs not described in the installation, operating and maintenance instructions must not be performed without consulting the manufacturer first.

⚠ DANGER

Strictly observe the safety datasheets or the safety regulations that are valid for the media used.

In cases of uncertainty:

- x Consult the nearest GEMÜ sales office.

2.2 Warning notes

Wherever possible, warning notes are organised according to the following scheme:

⚠ SIGNAL WORD

Type and source of the danger

- Possible consequences of non-observance.
- Measures for avoiding danger.

Warning notes are always marked with a signal word and sometimes also with a symbol for the specific danger. The following signal words and danger levels are used:

⚠ DANGER

Imminent danger!

- Non-observance will lead to death or severe injury.

⚠ WARNING

Potentially dangerous situation!

- Non-observance can cause death or severe injury.

⚠ CAUTION

Potentially dangerous situation!

- Non-observance can cause moderate to light injury.

CAUTION (WITHOUT SYMBOL)

Potentially dangerous situation!

- Non-observance can cause damage to property.

2.3 Symbols used

 Danger - hot surfaces!

 Danger - corrosive materials!

 Hand: indicates general information and recommendations.

●	Bullet point: indicates the tasks to be performed.
➤	Arrow: indicates the response(s) to tasks.
x	Enumeration sign

### 3 Definition of terms

#### Working medium

The medium that flows through the diaphragm valve.

#### Control medium

The medium whose increasing or decreasing pressure causes the diaphragm valve to be actuated and operated.

#### Control function

The possible actuation functions of the diaphragm valve.

### 4 Intended area of use

- x The GEMÜ 650 diaphragm valve is designed for installation in piping systems. It controls a flowing medium by being closed or opened by a control medium.

- x The valve may only be used providing the product technical criteria are complied with (see chapter 5 "Technical data").
- x Do not paint the bolts and plastic parts of the diaphragm valve!

#### ⚠ WARNING

**Use the diaphragm valve only for the intended purpose!**

- Otherwise the manufacturer liability and guarantee will be void.
- Use the diaphragm valve only in accordance with the operating conditions specified in the contract documentation and in the installation, operating and maintenance instructions.
- The diaphragm valve may only be used in potentially explosive zones confirmed in the declaration of conformity (ATEX).

## 5 Technical data

#### Working medium

Corrosive, inert, gaseous and liquid media which have no negative impact on the physical and chemical properties of the body and diaphragm material.

The valve will seal in both flow directions up to full operating pressure (gauge pressure).

#### Temperatures

**Medium temperature** -10 to 100 °C

##### Sterilisation temperature <sup>(1)</sup>

EPDM (code 13/3A)	max. 150 °C <sup>(2)</sup> , max. 60 min per cycle
EPDM (code 17)	max. 150 °C <sup>(2)</sup> , max. 180 min per cycle
EPDM (code 19)	max. 150 °C <sup>(2)</sup> , max. 180 min per cycle
EPDM (code 36)	max. 150 °C <sup>(2)</sup> , max. 60 min per cycle
PTFE/EPDM (code 54)	max. 150 °C <sup>(2)</sup> , no time limit per cycle
PTFE/EPDM (code 5M, 5Q)	max. 150 °C <sup>(2)</sup> , no time limit per cycle
PTFE/PVDF/EPDM (code 71)	not applicable

<sup>1</sup> The sterilisation temperature is valid for steam (saturated steam) or superheated water.


<sup>2</sup> If the sterilisation temperatures listed above are applied to the EPDM diaphragms for longer periods of time, the service life of the diaphragms will be reduced. In these cases, maintenance cycles must be adapted accordingly. This also applies to PTFE diaphragms exposed to high temperature fluctuations.

PTFE diaphragms can also be used as steam barriers; however, this will reduce their service life.

The maintenance cycles must be adapted accordingly.

GEMÜ 555 and 505 globe valves are particularly suitable for use in the area of steam generation and distribution.

The following valve arrangement for interfaces between steam pipes and process pipes has proven itself over time: A globe valve for shutting off steam pipes and a diaphragm valve as an interface to the process pipes.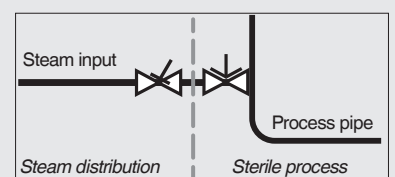

**Ambient temperature**

0 ... 60 °C

Control medium						
Inert gases						
Max. permissible temperature of control medium						60 °C
Filling volume						
Diaphragm size	DN	Actuator size	Actuator version	Spring set	Control function 1	Control function 2
8	4 to 15	0	T/R	1	0.01 dm <sup>3</sup>	0.01 dm <sup>3</sup>
			T/R	A	0.02 dm <sup>3</sup>	0.01 dm <sup>3</sup>
10	10 to 20	1	T/R/D/B	1	0.03 dm <sup>3</sup>	0.07 dm <sup>3</sup>
25	15 to 25	2	T/R/D/B	1	0.13 dm <sup>3</sup>	0.22 dm <sup>3</sup>
40	32 to 40	3	T/R/D/B	1	0.23 dm <sup>3</sup>	0.50 dm <sup>3</sup>
			T/R	A	0.50 dm <sup>3</sup>	-
50	50 to 65	4	T/R/D/B	1	0.50 dm <sup>3</sup>	1.20 dm <sup>3</sup>
80	65 to 80	5	T/R	1	2.68 dm <sup>3</sup>	3.20 dm <sup>3</sup>
			T/R	A/B	2.13 dm <sup>3</sup>	-
100	100	6	T/R	1	2.78 dm <sup>3</sup>	3.40 dm <sup>3</sup>
			T/R	A	2.15 dm <sup>3</sup>	-
150	150	8	T	A	5.30 dm <sup>3</sup>	-
C.f. 3 = for filling volume in open position see c.f. 1; for filling volume in closed position see c.f. 2						

Operating pressure [bar]								
MG	DN	Control function	Actuator version	EPDM		PTFE		
				Diaphragm material	All valve body materials	Diaphragm material	Forged and block material bodies	Investment cast body
8	4 to 15	1	0T1, 0R1	13, 17, 19, 36	0 to 8	54	0 to 6	0 to 6
			0TA, 0RA		0 to 10		0 to 10	0 to 6
		2 + 3	0T1, 0R1, 0TA, 0RA		0 to 10		0 to 10	0 to 6
							0 to 10	0 to 6
10	10 to 20	1	1T1, 1R1	13, 17, 19, 36	0 to 10	54	0 to 10	0 to 6
			1D1, 1B1		0 to 10		0 to 6	0 to 6
		2 + 3	1T1, 1R1		0 to 10		0 to 10	0 to 6
			1D1, 1B1		0 to 10		0 to 6	0 to 6
25	15 to 25	1	2T1, 2R1	13, 17, 19, 36	0 to 10	54, 5M	0 to 10	0 to 6
			2D1, 2B1		0 to 10		0 to 6	0 to 6
		2 + 3	2T1, 2R1		0 to 10		0 to 10	0 to 6
			2D1, 2B1		0 to 10		0 to 6	0 to 6
40	32 to 40	1	3T1, 3R1, 3D1, 3B1	13, 17, 19, 36	0 to 10	54, 5M	0 to 6	0 to 6
			3TA, 3RA		-		0 to 10	0 to 6
		2 + 3	3T1, 3R1		0 to 10		0 to 10	0 to 6
			3D1, 3B1		0 to 10		0 to 6	0 to 6
50	50 to 65	1	4T1, 4R1	13, 17, 19, 36	0 to 10	54, 5M	0 to 10	0 to 6
			4D1, 4B1		0 to 10		0 to 6	0 to 6
		2 + 3	4T1, 4R1		0 to 10		0 to 10	0 to 6
			4D1, 4B1		0 to 10		0 to 6	0 to 6
80	65 to 80	1	5T1, 5R1	13, 17, 19, 36	0 to 8	54, 5M	0 to 5	-
			5TA, 5RA		-		0 to 10	-
			5TB, 5RB		0 to 10		-	-
		2 + 3	5T1, 5R1		0 to 10		0 to 10	-
100	100	1	6T1, 6R1	13, 17, 19, 36	0 to 6	54, 5M	0 to 4	-
			6TA, 6RA		0 to 10		0 to 10	-
		2 + 3	6T1, 6R1		0 to 10		0 to 10	-
150	150	1	8TA, 8RA	-	-	5Q	0 to 10	-

All pressures are gauge pressures. Operating pressure values were determined with static operating pressure applied on one side of a closed valve. Sealing at the valve seat and atmospheric sealing is ensured for the given values.  
Information on operating pressures applied on both sides and for high purity media on request  
MG = diaphragm size

Control pressure [bar]				
MG	DN	Control function	Actuator version	Control pressure
8	4 to 15	1	0T1, 0R1	5.0 to 7.0
			0TA, 0RA	3.5 to 7.0
		2 + 3	0T1, 0R1	max. 5.5
			0TA, 0RA	max. 4.5
10	10 to 20	1	1T1, 1R1, 1D1, 1B1	4.5 to 7.0
		2 + 3	1T1, 1R1, 1D1, 1B1	max. 4.5
25	15 to 25	1	2T1, 2R1, 2D1, 2B1	5.0 to 7.0
		2 + 3	2T1, 2R1, 2D1, 2B1	max. 4.5
40	32 to 40	1	3T1, 3R1, 3D1, 3B1	4.5 to 7.0
			3TA, 3RA	3.5 to 7.0
		2 + 3	3T1, 3R1, 3D1, 3B1	max. 4.5
50	50 to 65	1	4T1, 4R1, 4D1, 4B1	4.5 to 7.0
		2 + 3	4T1, 4R1, 4D1, 4B1	max. 4.5
80	65 to 80	1	5T1, 5R1	3.5 to 7.0
			5TA, 5RA	4.5 to 7.0
			5TB, 5RB	4.0 to 7.0
		2 + 3	5T1, 5R1	max. 4.0
100	100	1	6T1, 6R1	3.5 to 7.0
			6TA, 6RA	5.0 to 7.0
		2 + 3	6T1, 6R1	max. 4.0
150	150	1	8TA, 8RA	7.0 to 8.0

MG = diaphragm size

Kv values [m³/h]									
Pipe standard	DIN	EN 10357 series B (formerly DIN 11850 series 1)	EN 10357 series A (formerly DIN 11850 series 2) / DIN 11866 series A	DIN 11850 Series 3	SMS 3008	ASME BPE / DIN 11866 series C	ISO 1127 / EN 10357 series C / DIN 11866 series B	DIN ISO 228	
Connection code	0	16	17	18	37	59	60	1	
MG	DN								
8	4	0.5	-	-	-	-	-	-	-
	6	-	-	1.1	-	-	-	1.2	-
	8	-	-	1.3	-	-	0.6	2.2	1.4
	10	-	2.1	2.1	2.1	-	1.3	-	-
	15	-	-	-	-	-	2.0	-	-
10	10	-	2.4	2.4	2.4	-	2.2	3.3	-
	12	-	-	-	-	-	-	-	3.2
	15	3.3	3.8	3.8	3.8	-	2.2	4.0	3.4
	20	-	-	-	-	-	3.8	-	-
25	15	4.1	4.7	4.7	4.7	-	-	7.4	6.5
	20	6.3	7.0	7.0	7.0	-	4.4	13.2	10.0
	25	13.9	15.0	15.0	15.0	12.6	12.2	16.2	14.0
40	32	25.3	27.0	27.0	27.0	26.2	-	30.0	26.0
	40	29.3	30.9	30.9	30.9	30.2	29.5	32.8	33.0
50	50	46.5	48.4	48.4	48.4	51.7	50.6	55.2	60.0
	65	-	-	-	-	62.2	61.8	-	-
80	65	-	-	77.0	-	68.5	68.5	96.0	-
	80	-	-	111.0	-	80.0	87.0	111.0	-
100	100	-	-	194.0	-	173.0	188.0	214.0	-
150	150	-	-	-	-	-	570.0	-	-

MG = diaphragm size

Kv values determined acc. to DIN EN 60534, inlet pressure 5 bar,  $\Delta p$  1 bar, stainless steel valve body (forged body) and soft elastomer diaphragm.


The Kv values for other product configurations (e.g. other diaphragm or body materials) may differ. In general, all diaphragms are subject to the influences of pressure, temperature, the process and their tightening torques. Therefore the Kv values may exceed the tolerance limits of the standard.

The Kv value curve (Kv value dependent on valve stroke) can vary depending on the diaphragm material and duration of use.


## Autoclavability

Actuator size 0	Standard version with autoclave capability
Actuator size 1	Standard version with autoclave capability
Actuator size 2	Standard version with autoclave capability
Actuator size 3	with special version
Actuator size 4	with special version
Actuator size 5	not possible
Actuator size 6	not possible
Actuator size 8	not possible


Control function 2 + 3  
with elastomer diaphragm  
Diaphragm size 8 to 25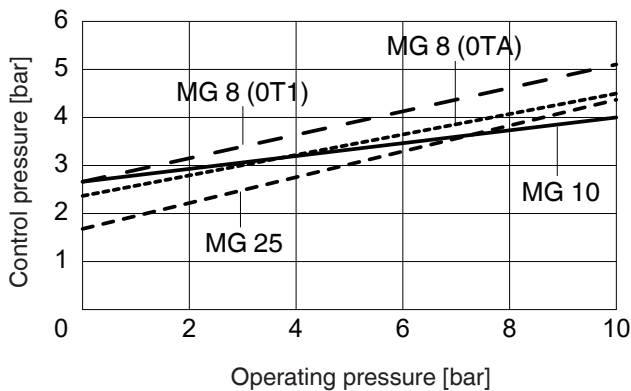


Control function 2 + 3  
with elastomer diaphragm  
Diaphragm size 40 to 100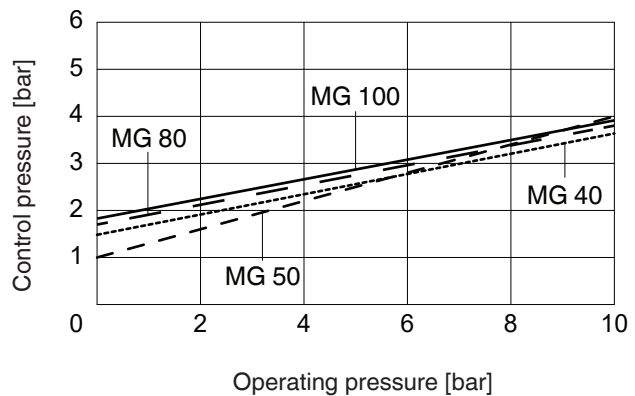

Control function 2 + 3  
with PTFE diaphragm  
Diaphragm size 8 to 25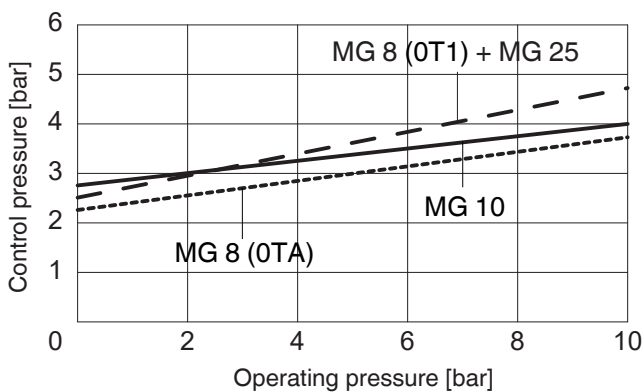

Control function 2 + 3  
with PTFE diaphragm  
Diaphragm size 40 to 100

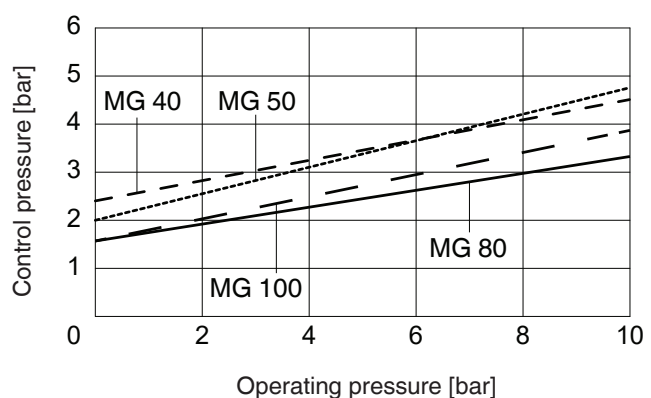The control pressure depending on the prevailing operating pressure, as shown in the diagram, is intended as a guide for operating the system with low wear on the diaphragm.

## 6 Order data (2/2-way valves)

Body configuration	Code
Tank bottom valve body (actuator version T)	B**
2/2-way body (actuator version D and T)	D
T body (actuator version T)	T*
* For dimensions see T Valves brochure	
** Dimensions and versions on request	

Connection	Code
<b>Butt weld spigots</b>	
Spigots DIN	0
Spigots EN 10357 series B (formerly DIN 11850 series 1)	16
Spigot EN 10357 series A (formerly DIN 11850 series 2) / DIN 11866 series A	17
Spigots DIN 11850 series 3	18
Spigots JIS-G 3447	35
Spigots JIS-G 3459	36
Spigots SMS 3008	37
Spigots BS 4825 Part 1	55
Spigot ASME BPE / DIN 11866 series C	59
Spigot ISO 1127 / EN 10357 series C / DIN 11866 series B	60
Spigots ANSI/ASME B36.19M Schedule 10s	63
Spigots ANSI/ASME B36.19M Schedule 5s	64
Spigots ANSI/ASME B36.19M Schedule 40s	65
<b>Threaded connections</b>	
Threaded sockets DIN ISO 228	1
Threaded spigots DIN 11851	6
Cone spigot and union nut DIN 11851	6K
Aseptic unions on request	
<b>Flanges</b>	
Flanges EN 1092 / PN16 / form B, length EN 558, series 1, ISO 5752, basic series 1	8*
Flanges ANSI Class 150 RF, length MSS SP-88	38*
Flanges ANSI Class 125/150 RF, length EN 558, series 1, ISO 5752, basic series 1	39*
<b>Clamp connections</b>	
Clamps ASME BPE for pipe ASME BPE, length ASME BPE	80
Clamps DIN 32676 series B for pipe EN ISO 1127, length EN 558, series 7	82
Clamps ASME BPE for pipe ASME BPE, length EN 558, series 7	88
Clamps DIN 32676 series A for pipe DIN 11850, length EN 558, series 7	8A
Clamps SMS 3017 for pipe SMS 3008, length EN 558, series 7	8E
Clamps DIN 32676 series C, length FTF ASME BPE	8P
Clamps DIN 32676 series C, length FTF EN 558 series 7	8T
Aseptic clamps on request	
* Connection code 8, 38, 39 only possible in conjunction with actuator version code B / R	

Valve body material	Code
1.4435, investment casting	C3
1.4408, investment casting	37
1.4408, PFA lined	39
1.4435 (316L), forged body	40
1.4435 (316L), block material	41
1.4435 (BN2), forged body $\Delta$ Fe<0.5%	42
1.4435 (BN2), block material, $\Delta$ Fe<0,5 %	43
1.4539, forged body	F4

Diaphragm material	Code
EPDM	13 3A*
EPDM	17
EPDM	19
EPDM	36
PTFE/EPDM, one-piece	54
PTFE/EPDM, two-piece	5M**
PTFE/EPDM, two-piece	5Q
PTFE/PVDF/EPDM, three-piece	71***
* for diaphragm size 8	
** Code 5M not available in diaphragm size 10	
*** Code 71 only available for bodies with PFA lining (Code 39)	
Material complies with FDA requirements	

Control function	Code
Normally closed (NC)	1
Normally open (NO)	2
Double acting (DA) (with opening spring)	3

Actuator size	Code
Actuator size 0 (diaphragm size 8)	0
Actuator size 1 (diaphragm size 10)	1
Actuator size 2 (diaphragm size 25)	2
Actuator size 3 (diaphragm size 40)	3
Actuator size 4 (diaphragm size 50)	4
Actuator size 5 (diaphragm size 80)	5
Actuator size 6 (diaphragm size 100)	6
Actuator size 8 (diaphragm size 150)	8

Actuator version	Code
Only for body configuration D (diaphragm size 10 - 50)	D
For body config. D (diaphragm size 10 - 50) Control air connector 90° offset to flow direction	B
For body config. B, D, M and T (diaphragm size 8 - 100)	T
For body config. B, D, M and T (diaphragm size 8 - 100) Control air connector 90° offset to flow direction	R

Spring set	Code
Standard	1
For higher operating pressure	A
For higher operating pressure	B

## Internal surface finishes for forged and block material bodies <sup>1)</sup>

Readings for Process Contact Surfaces	Mechanically polished <sup>2)</sup>		Electropolished	
	Hygienic class DIN 11866	Code	Hygienic class DIN 11866	Code
Ra ≤ 0.80 µm	H3	1502	HE3	1503
Ra ≤ 0.60 µm	-	1507	-	1508
Ra ≤ 0.40 µm	H4	1536	HE4	1537
Ra ≤ 0.25 µm <sup>3)</sup>	H5	1527	HE5	1516

Readings for Process Contact Surfaces acc. to ASME BPE 2016 <sup>4)</sup>	Mechanically polished <sup>2)</sup>		Electropolished	
	ASME BPE Surface Designation	Code	ASME BPE Surface Designation	Code
Ra Max. = 0.76 µm (30 µinch)	SF3	SF3	-	-
Ra Max. = 0.64 µm (25 µinch)	SF2	SF2	SF6	SF6
Ra Max. = 0.51 µm (20 µinch)	SF1	SF1	SF5	SF5
Ra Max. = 0.38 µm (15 µinch)	-	-	SF4	SF4

## Internal surface finishes for investment cast bodies

Readings for Process Contact Surfaces	Mechanically polished <sup>2)</sup>	
	Hygienic class DIN 11866	Code
Ra ≤ 6.30 µm	-	1500
Ra ≤ 0.80 µm	H3	1502
Ra ≤ 0.60 µm <sup>5)</sup>	-	1507

<sup>1)</sup> Surface finishes of customized valve bodies may be limited in special cases.

<sup>2)</sup> Or any other finishing method that meets the Ra value (acc. to ASME BPE).

<sup>3)</sup> The smallest possible Ra finish for pipe connections with an internal pipe diameter < 6 mm is 0.38 µm.

<sup>4)</sup> When using these surfaces, the bodies are marked according to the specifications of ASME BPE.

The surfaces are only available for valve bodies which are made of materials (e.g. GEMÜ material codes 40, 41, F4, 44) and use connections (e.g. GEMÜ connection codes 59, 80, 88) according to ASME BPE.

<sup>5)</sup> Not possible for GEMÜ connection code 59, DN 8 and GEMÜ connection code 0, DN 4.

Ra acc. to DIN EN ISO 4288 and ASME B46.1

Special function	Code
3-A compliant design	M

Order example	650	50	D	60	40	54	1	4	T	1	1503	M
Type	650											
Nominal size		50										
Body configuration (code)			D									
Connection (code)				60								
Valve body material (code)					40							
Diaphragm material (code)						54						
Control function (code)							1					
Actuator size (code)								4				
Actuator version (code)									T			
Spring set (code)										1		
Surface finish (code)											1503	
Special function (Code)												M

## 7 Manufacturer's information

### 7.1 Transport

- Only transport the diaphragm valve by suitable means. Do not drop. Handle carefully.
- Dispose of packaging material according to relevant local or national disposal regulations/environmental protection laws.

### 7.2 Delivery and performance

- Check that all parts are present and check for any damage immediately upon receipt.
- The scope of delivery is apparent from the dispatch documents and the design from the order number.
- Condition of the valve as supplied to customer:

Control function:	Condition:
1 Normally closed (NC)	closed
2 Normally open (NO)	open
3 Double acting (DA)	open

- The performance of the diaphragm valve is checked at the factory.

### 7.3 Storage

- Store the diaphragm valve free from dust and moisture in its original packaging.
- Avoid UV rays and direct sunlight.
- Maximum storage temperature: 40 °C.
- Solvents, chemicals, acids, fuels or similar fluids must not be stored in the same room as valves and their spare parts.


### 7.4 Tools required

- The tools required for installation and assembly are **not** included in the scope of delivery.
- Use appropriate, functional and safe tools.

## 8 Functional description

GEMÜ 650 is a metal diaphragm valve with 2/2-way, T or tank bottom valve body or multi-port design. The valve has a piston actuator and an optical position indicator as standard. All metallic actuator components are made of stainless steel. The compression springs of diaphragm sizes 80 and 100 are made of epoxy coated spring steel. "Normally closed (NC)", "Normally open (NO)" and "Double acting (DA)" control functions are available. The valve body and the diaphragm are available in various designs as shown in the datasheet. The valve can be cleaned (CIP) and sterilized (SIP) without dismantling (autoclavable depending on version). Diverse accessories are available, e.g. stroke limiters, electrical position indicators, positioners.


## 9 Construction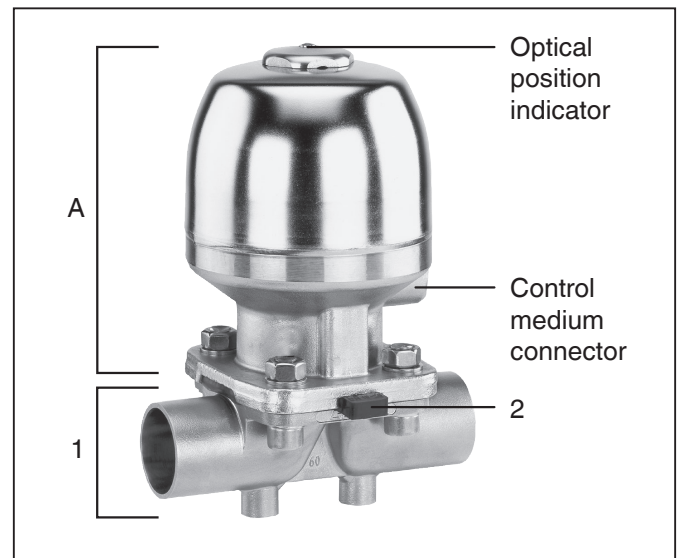

Construction

1	Valve body
2	Diaphragm
A	Actuator

## 9.1 Product label

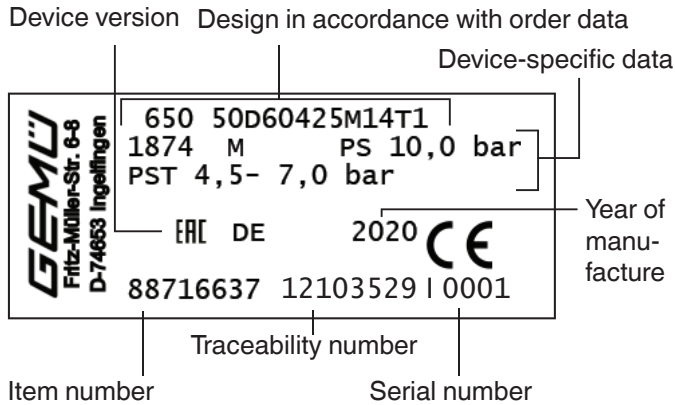The month of manufacture is encoded in the traceability number and can be obtained from GEMÜ.

The product was manufactured in Germany.

## 10 Installation and operation

### Prior to installation:

- Ensure that valve body and diaphragm material are appropriate and compatible to handle the working medium.
- **Check the suitability prior to the installation.**  
See chapter 5 "Technical data".

### 10.1 Installing the diaphragm valve

#### ⚠ WARNING

##### The equipment is subject to pressure!

- Risk of severe injury or death!
- Only work on depressurized plant.

#### ⚠ WARNING

##### Corrosive chemicals!

- Risk of caustic burns!
- Wear appropriate protective gear when installing.

#### ⚠ CAUTION

##### Hot plant components!

- Risk of burns!
- Only work on plant that has cooled down.

#### ⚠ CAUTION

##### Never use the valve as a step or an aid for climbing!

- This entails the risk of slipping-off or damaging the valve.

#### CAUTION

##### Do not exceed the maximum permissible pressure!

- Take precautionary measures to avoid possible pressure surges (water hammer).

- Installation work must only be performed by trained personnel.
- Use appropriate protective gear as specified in plant operator's guidelines.

### Installation location:

#### ⚠ CAUTION

- Do not apply external force to the valve.
- Choose the installation location so that the valve cannot be used as a foothold.
- Lay the pipeline so that the valve body is protected against transverse and bending forces, and also vibrations and tension.
- Only install the valve between matching aligned pipes.

- x Direction of the working medium: optional.
- x Mounting position of the diaphragm valve: optional.

### Installation:

1. Ensure the suitability of the valve for each respective use. The valve must be appropriate for the piping system operating conditions (medium, medium concentration, temperature and pressure) and the prevailing ambient conditions. Check the technical data of the valve and the materials.
2. Shut off plant or plant component.
3. Secure against recommissioning.
4. Depressurize the plant or plant component.

5. Completely drain the plant or plant component and allow it to cool down until the temperature is below the media vaporization temperature and cannot cause scalding.
6. Correctly decontaminate, rinse and ventilate the plant or plant component.

#### **Note on EHEDG certified valves:**

EHEDG certified valves must be installed so that they can be easily cleaned and drained. For valves with weld ends, the weld seams must be made in accordance with EHEDG Guideline 9 and 35. For valves with removable connections, the "Position Paper" of the EHEDG must be taken into account and, if necessary, special seals must be used.

#### **Installation - Butt weld spigots:**

1. Adhere to good welding practices!
2. Disassemble the actuator with the diaphragm before welding the valve body into the pipeline (see chapter 11.1).
3. Allow the butt weld spigots to cool down.
4. Reassemble the valve body and the actuator with diaphragm (see chapter 11.4).

#### **Installation - Clamp connections:**

- When installing clamp connections, insert a gasket between the body clamp and the adjacent piping clamp and join them using the appropriate clamp fitting. The gasket and the clamp for clamp connections are not included in the scope of delivery.

#### **Important:**


Butt weld spigots / clamp connections: Angle of rotation for welding into pipeline to enable optimised draining see brochure "Angle of rotation for 2/2-way valve bodies" (on request or [www.gemu-group.com](http://www.gemu-group.com)).

#### **Installation - Threaded connections:**

- Screw the threaded connections into the piping in accordance with valid standards.
- Screw the diaphragm valve body into the piping, use appropriate thread sealant. The thread sealant is not included in the scope of delivery.

#### **Installation - Flanges:**

1. Pay attention to clean, undamaged sealing surfaces on the mating flanges.
2. Align flanges carefully before installing them.
3. Centre the seals accurately.
4. Connect the valve flange and the piping flange using appropriate sealing material and matching bolting. Sealing material and bolts are not included in the scope of delivery.
5. Use all flange holes.
6. Only use connector elements made of approved materials!
7. Tighten the bolts diagonally!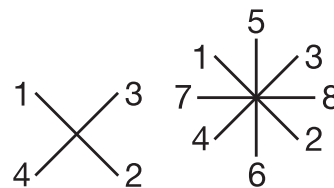

#### **Observe appropriate regulations for connections!**

#### **After the installation:**

- Re-attach or reactivate all safety and protective devices.

## **10.2 Control functions**

The following control functions are available:

#### **Control function 1**

##### **Normally closed (NC):**

Valve resting position: closed by spring force. Activation of the actuator (connector 2) opens the valve. When the actuator is vented, the valve is closed by spring force.

#### **Control function 2**


##### **Normally open (NO):**

Valve resting position: opened by spring force. Activation of the actuator (connector 4) closes the valve. When the actuator is vented, the valve is opened by spring force.


**Control function 3**  
**Double acting (DA):**

Valve resting position: opened by spring force. The valve is opened and closed by activating the respective control medium connectors (connector 2: open / connector 4: close).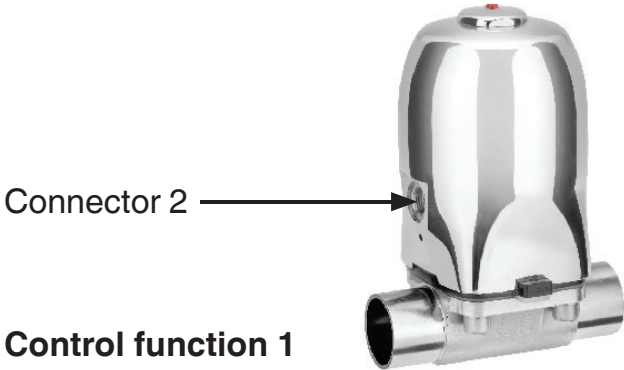

**Control function 1**

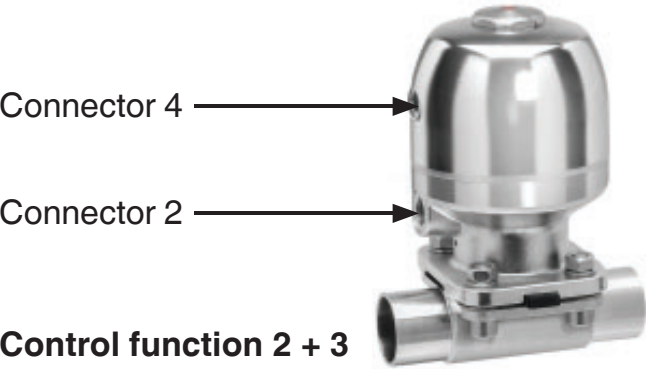**Control function 2 + 3**


Thread size of the control medium connectors:

Diaphragm size 8: G1/8

Diaphragm size 10 - 100: G1/4

Control function		Connectors
1	Normally closed (NC)	2: Control medium (open)
2	Normally open (NO)	4: Control medium (close)
3	Double acting (DA)	2: Control medium (open) 4: Control medium (close)
For connectors 2 / 4 see pictures on page 11		

**10.4 Optical position indicator**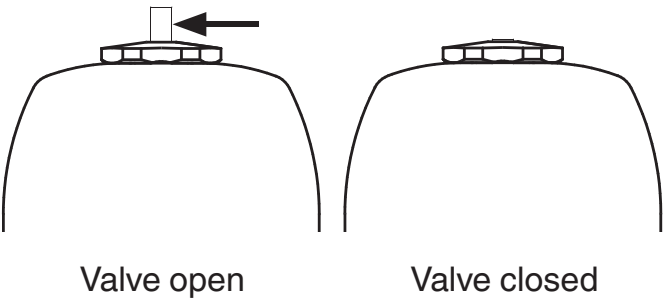

Control function	Connectors	
	2	4
1 (NC)	+	-
2 (NO)	-	+
3 (DA)	+	+
+ = available / - = not available (for connectors 2 / 4 see pictures above)		


**10.3 Connecting the control medium**

**Important:**

Connect the control medium lines tension-free and without any bends or knots!


Use appropriate connectors according to the application.

## 11 Fitting/removing spare parts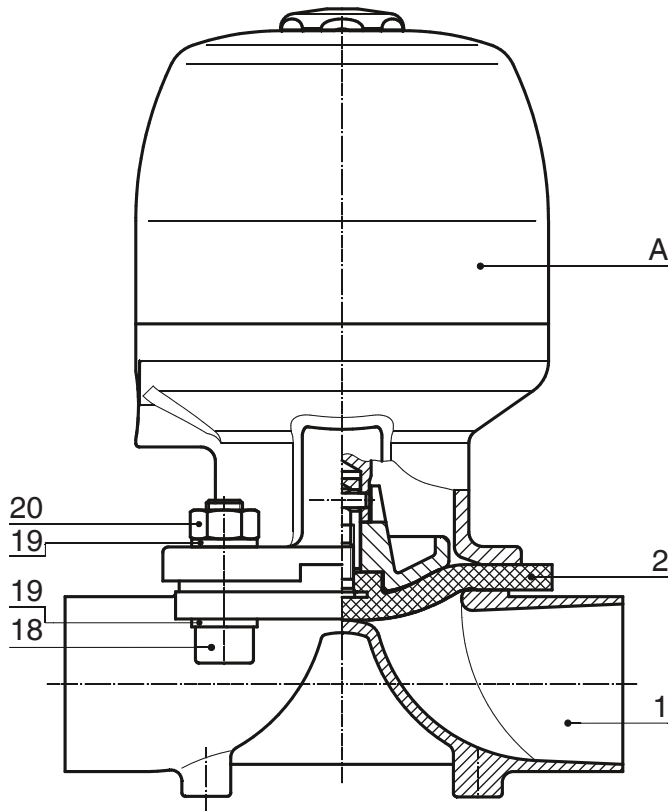

### 11.1 Valve disassembly (removing actuator from body)

1. Move actuator **A** to the open position.
2. Loosen the fastening elements between the valve body **1** and the actuator **A** diagonally and remove them.

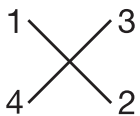3. Lift actuator **A** off valve body **1**.
4. Move actuator **A** to the closed position.

#### **Important:**

After disassembly, clean all parts of contamination (do not damage parts). Check parts for potential damage, replace if necessary (only use genuine parts from GEMÜ).

## 11.2 Removing the diaphragm

#### **Important:**

Before removing the diaphragm, please remove the actuator, see "Valve disassembly (removing actuator from body)".

1. Unscrew the diaphragm or pull it out (diaphragm size 8).
2. Clean all parts of remains of product and contaminants. Do not scratch or damage parts during cleaning!
3. Check all parts for potential damage.
4. Replace damaged parts (only use genuine parts from GEMÜ).

## 11.3 Mounting the diaphragm

### 11.3.1 General information

#### **Important:**

Mount the correct diaphragm that suits the valve (suitable for medium, medium concentration, temperature and pressure). The diaphragm is a wearing part. Check the technical condition and function of the diaphragm valve before commissioning and during the whole duration of use. Carry out checks regularly and determine the check intervals in accordance with the conditions of use and/or the regulatory codes and provisions applicable for this application.

**Important:**

If the diaphragm is not screwed into the adapter far enough, the closing force is transmitted directly onto the diaphragm pin and not via the compressor. This will cause damage and early failure of the diaphragm and thus leakage of the valve. If the diaphragm is screwed in too far, perfect sealing at the valve seat will not be achieved. The function of the valve is no longer ensured.

**Important:**

Incorrectly mounted diaphragm may cause valve leakage / emission of medium. In this case remove the diaphragm, check the complete valve and diaphragm and reassemble again proceeding as described above.

Diaphragm sizes 10 - 80:


The compressor is loose.

Diaphragm sizes 8 and 100:

The compressor is fixed to the spindle.


**Diaphragm size 8:**

Compressor and actuator flange seen from below:

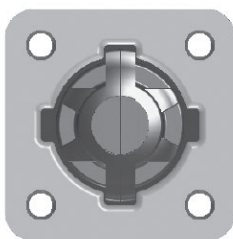
**Diaphragm size 10:**

Compressor and actuator flange seen from below:

Picture 1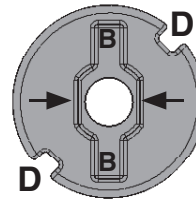

Picture 2

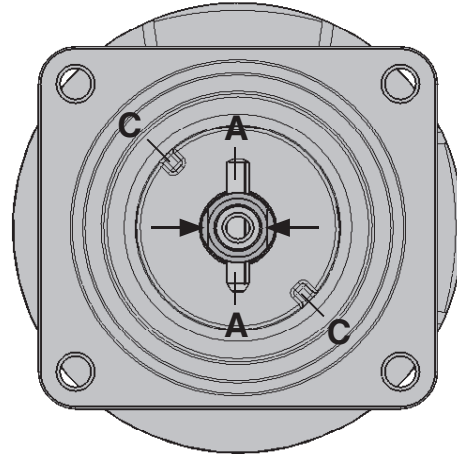
**Anti-twist system of the spindle at the compressor**

A double flat (arrows picture 2) is fitted at the end of the actuator spindle to protect the actuator spindle against twisting. When mounting the compressor, the double flat must be in correct alignment with the recess of the compressor back (arrows picture 1). If the actuator spindle is not in the correct position, it must be turned to the correct position. The position of **A** is offset by 45° to the position of **C**.

Place the compressor loosely on the actuator spindle, fit the recesses **D** into the guides **C** and **A** into **B**. The compressor must be able to be moved freely between the guides!

### Diaphragm size 25 - 80:


Compressor and actuator flange seen from below: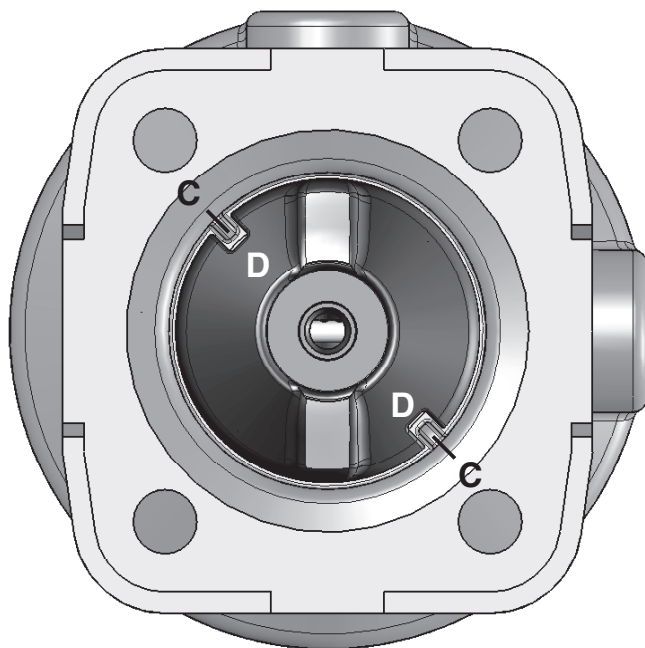

Place the compressor loosely on the actuator spindle, fit the recesses **D** into the guides **C**. The compressor must be able to be moved freely between the guides!

### Diaphragm size 100:


Compressor and actuator flange seen from below: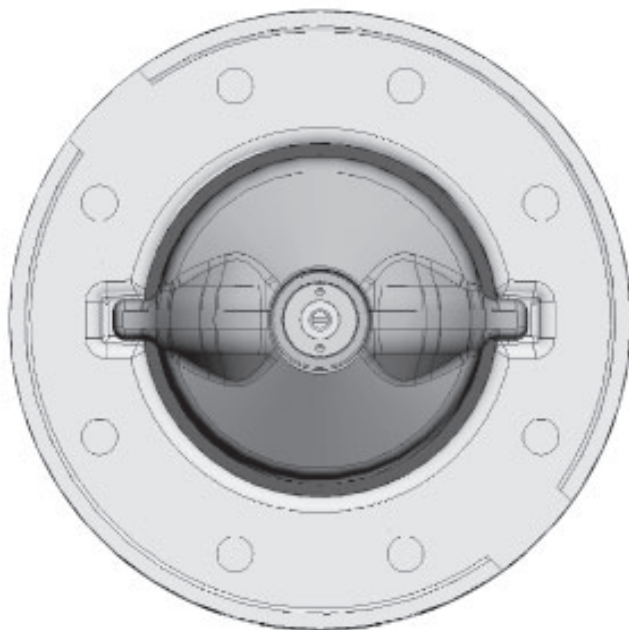

## 11.3.2 Mounting a concave diaphragm

### Diaphragm size 8

#### Push-fit diaphragm:

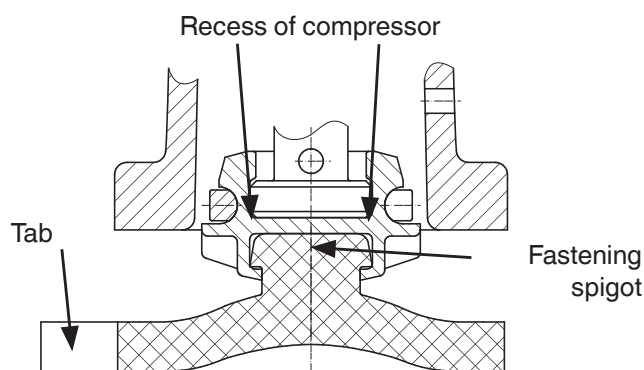1. Move actuator **A** to the closed position.
2. Place the diaphragm **2** with the fastening spigot in an inclined position at the recess of the compressor.


#### Important:

Do not use greases or lubricants!

3. Turn the diaphragm as manual force is applied to push the spigot into the compressor.
4. Align diaphragm tab with identifying manufacturer and material in parallel to compressor weir.

### Diaphragm sizes 10 - 100

#### Threaded pin type diaphragm: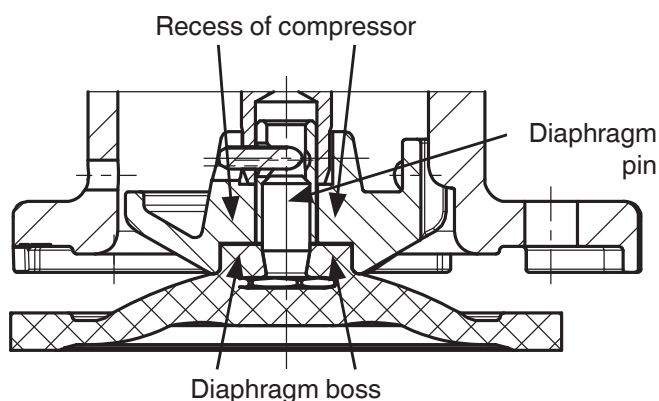


1. Move actuator **A** to the closed position.
2. Diaphragm size 10: Ensure that the anti-twist system is engaged.  
Diaphragm sizes 25 - 80: Place the compressor loosely on the actuator spindle, fit the recesses into the guides (see chapter 11.3.1 "General information")


3. Check if the compressor fits closely in the guides.
4. Screw new diaphragm tightly into the compressor manually.
5. Check if the diaphragm boss fits closely in the recess of the compressor.
6. If it is difficult to screw it in, check the thread, replace damaged parts (only use genuine parts from GEMÜ).
7. When clear resistance is felt turn back the diaphragm anticlockwise until its bolt holes are in correct alignment with the bolt holes of the actuator.

### 11.3.3 Mounting a convex diaphragm

1. Move actuator **A** to the closed position.
2. Diaphragm sizes 25 - 80: Place the compressor loosely on the actuator spindle, fit the recesses into the guides (see chapter 11.3.1 "General information")
3. Check if the compressor fits closely in the guides.
4. Invert the new diaphragm face manually; use a clean, padded mat with bigger nominal sizes.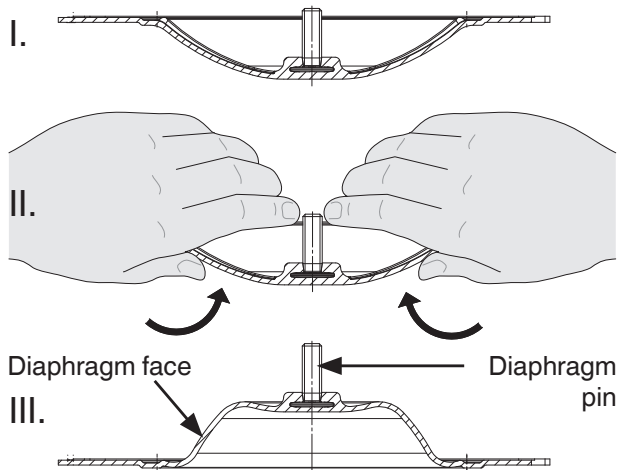


5. Position the new backing diaphragm onto the compressor.
6. Position the diaphragm face onto the backing diaphragm.
7. Screw diaphragm face tightly into the compressor manually. The diaphragm boss must fit closely in the recess of the compressor.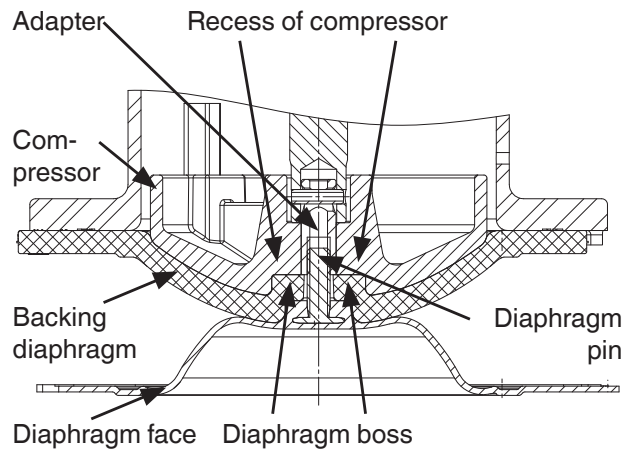

8. If it is difficult to screw it in, check the thread, replace damaged parts.
9. When clear resistance is felt turn back the diaphragm anticlockwise until its bolt holes are in correct alignment with the bolt holes of the actuator.
10. Press the diaphragm face tightly onto the backing diaphragm manually so that it returns to its original shape and fits closely on the backing diaphragm.

### 11.4 Actuator mounting on the valve body

1. Move actuator **A** to the open position.
2. Position actuator **A** with the mounted diaphragm **2** on the valve body **1**, take care to align the compressor weir and valve body weir (only for diaphragm size 8).
3. Tighten bolts **18**, washers **19** and nuts **20** by hand (hand tight only) (fastening elements may vary dependent on diaphragm size and / or valve body version).
4. Move actuator **A** to the closed position.
5. Fully tighten the bolts **18** with nuts **20** diagonally.

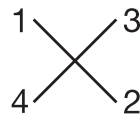6. Ensure that the diaphragm **2** is compressed evenly (approx. 10-15 %).  
⇒ Even compression is detected by an even bulge to the outside.

**Please note:** For a code 5M diaphragm (convex diaphragm), the PTFE diaphragm face and the EPDM backing diaphragm must be positioned level with and parallel to the valve body.

7. Check tightness of completely assembled valve.

	<p><b>Important:</b></p> <p>Service and maintenance: Diaphragms set in the course of time. After valve disassembly / assembly check that the bolts <b>18</b> and nuts <b>20</b> on the body are tight and retighten as necessary (at the very latest after the first sterilization process).</p>
---	--

## 12 Commissioning

⚠ WARNING	
	<p><b>Corrosive chemicals!</b></p> <ul style="list-style-type: none"> <li>➤ Risk of caustic burns!</li> <li>● Check the tightness of the media connections prior to commissioning!</li> <li>● Use only the appropriate protective gear when performing the tightness check.</li> </ul>

⚠ CAUTION
<p><b>Protect against leakage!</b></p> <ul style="list-style-type: none"> <li>● Provide precautionary measures against exceeding the maximum permitted pressures caused by pressure surges (water hammer).</li> </ul>

- Check the tightness and the function of the diaphragm valve (close and reopen the diaphragm valve).
- If the plant is new and after repairs rinse the piping system with a fully opened diaphragm valve (to remove any harmful foreign matter).

## 13 Inspection and maintenance

⚠ WARNING	
<p><b>The equipment is subject to pressure!</b></p> <ul style="list-style-type: none"> <li>➤ Risk of severe injury or death!</li> <li>● Only work on depressurized plant.</li> </ul>	

⚠ CAUTION	
	<p><b>Hot plant components!</b></p> <ul style="list-style-type: none"> <li>➤ Risk of burns!</li> <li>● Only work on plant that has cooled down.</li> </ul>

⚠ CAUTION	
<ul style="list-style-type: none"> <li>● Servicing and maintenance work must only be performed by trained personnel.</li> <li>● GEMÜ shall assume no liability whatsoever for damages caused by improper handling or third-party actions.</li> <li>● In case of doubt, contact GEMÜ prior to commissioning.</li> </ul>	

1. Use appropriate protective gear as specified in plant operator's guidelines.
  2. Shut off plant or plant component.
  3. Secure against recommissioning.
  4. Depressurize the plant or plant component.
- The operator must carry out regular visual examination of the valves dependent on the operating conditions and the potential danger in order to prevent leakage and damage. The valve also has to be disassembled in corresponding intervals and checked for wear (see chapter 11 "Fitting/removing spare parts").

## 14 Cleaning and sterilization

The valve can be cleaned (CIP) and sterilized (SIP) without removal. **The conditions in chapter "Technical data" (operating, cleaning and sterilization media, temperatures) must be observed.** During cleaning and sterilization, the valve must be permanently open.

## 15 Disassembly

Disassembly is performed observing the same precautionary measures as for installation.

- Disassemble the diaphragm valve (see chapter 11.1 "Valve disassembly (removing actuator from body)").

## 16 Disposal

- All valve parts must be disposed of according to relevant local or national disposal regulations / environmental protection laws.
- Pay attention to adhered residual material and gas diffusion from penetrated media.

## 17 Returns

- Clean the diaphragm valve.
- Request a return delivery note from GEMÜ.
- Returns must be made with a completed return delivery note.

If not completed, GEMÜ cannot process  
x credits or  
x repair work  
but will dispose of the goods at the operator's expense.

### Note for returns:

Legal regulations for the protection of the environment and personnel require that the completed and signed return delivery note is included with the dispatch documents. Returned goods can be processed only when this note is completed.

## 18 Information

### Note on Directive

#### 2014/34/EU (ATEX Directive):

A supplement to Directive 2014/34/EU is included with the product if it was ordered according to ATEX.

### Note on staff training:

Please contact us at the address on the last page for staff training information.

Should there be any doubts or misunderstandings in the preceding text, the German version of this document is the authoritative document!


## 19 Troubleshooting / Fault clearance

Error	Possible cause	Troubleshooting
Control medium escapes from vent hole*/vent* in the actuator cover for control function NC or connector 2 for control function NO (see chapter 10.2 "Control functions").	Piston faulty	Replace the actuator
Control medium escapes from leak detection hole*	Spindle seal leaking	Replace the actuator and check control medium for impurities
Working medium escapes from leak detection hole*	Shut off diaphragm faulty	Check shut off diaphragm for potential damage, replace diaphragm if necessary
Valve doesn't open or doesn't open fully	Control pressure too low (for control function NC)	Operate valve with control pressure specified in data sheet
	Pilot valve faulty	Check and replace pilot valve
	Actuator defective	Replace the actuator
	Control medium not connected	Connect the control medium
	Shut off diaphragm incorrectly mounted	Remove actuator, check diaphragm mounting, replace if necessary
	Actuator spring faulty (for control function NO)	Replace the actuator
Valve leaks downstream (doesn't close or doesn't close fully)	Operating pressure too high	Operate valve with operating pressure specified in datasheet
	Control pressure too low (for control function NO and control function DA)	Operate valve with control pressure specified in data sheet
	Foreign matter between shut off diaphragm and valve body weir	Remove actuator, remove foreign matter, check diaphragm and valve body weir for potential damage and replace if necessary
	Valve body weir leaking or damaged	Check valve body weir for damage, if necessary replace valve body
	Shut off diaphragm faulty	Check shut off diaphragm for potential damage, replace diaphragm if necessary
	Actuator spring faulty (for control function NC)	Replace the actuator
Valve leaks between actuator and valve body	Shut off diaphragm incorrectly mounted	Remove actuator, check diaphragm mounting, replace if necessary
	Bolting between valve body and actuator loose	Retighten bolting between valve body and actuator
	Shut off diaphragm faulty	Check shut off diaphragm for potential damage, replace diaphragm if necessary
	Actuator/valve body damaged	Replace the actuator/valve body
Valve body connection to piping leaks	Incorrect installation	Check installation of valve body in piping
	Unions/threaded connections loose	Tighten unions/threaded connections
	Sealing material faulty	Replace sealing material
Valve body leaks	Valve body faulty or corroded	Check valve body for potential damage, replace valve body if necessary

\* see chapter 19 "Sectional drawings and spare parts"


20 Sectional drawings and spare parts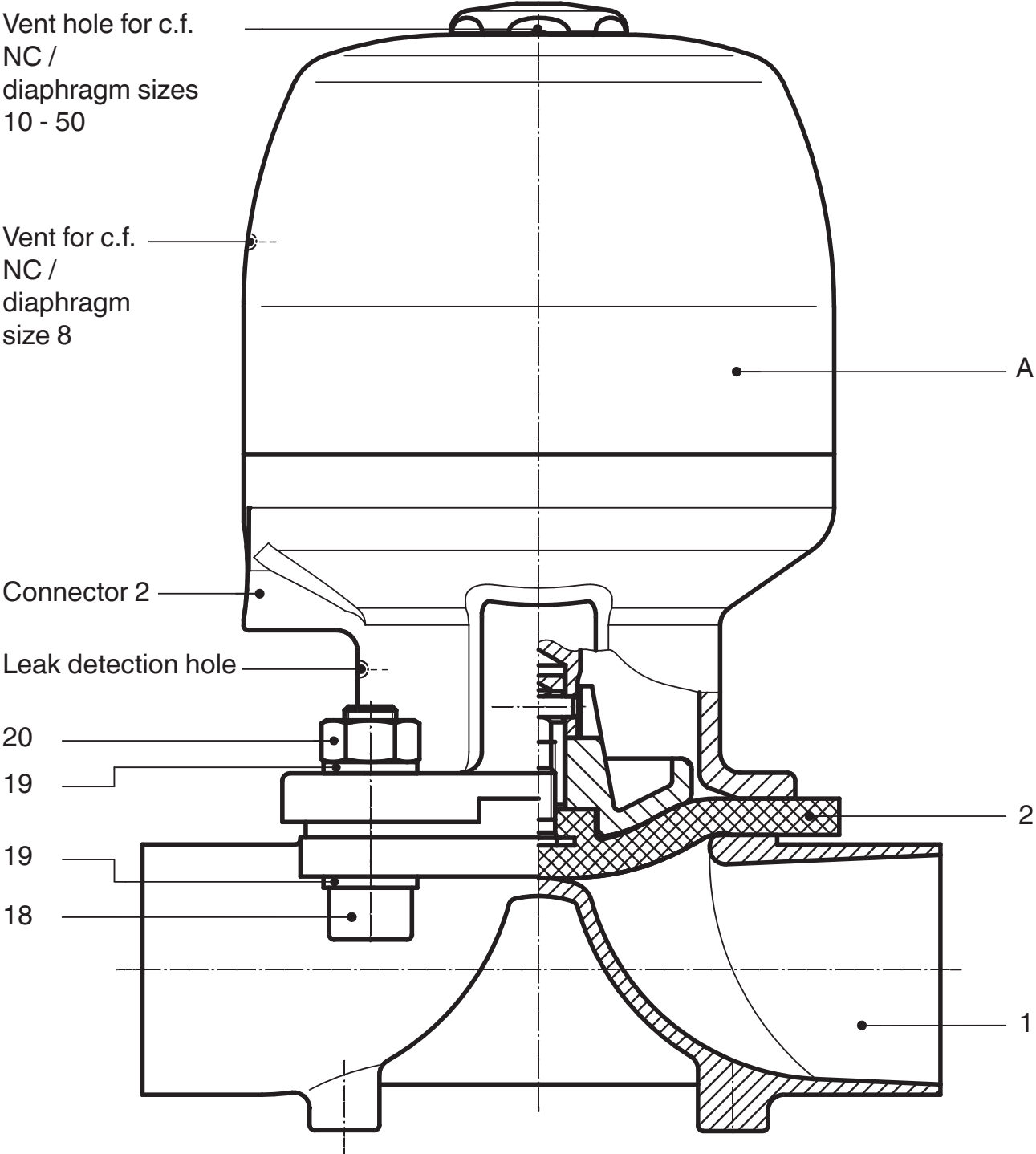


Item	Name	Order designation
Diaphragm sizes 8 - 50 / Actuator version: T		
1	Valve body	K600...
2	Diaphragm	600...M
18	Bolt	} 650...S30...
19	Washer	
20	Nut	
A	Actuator	9650...


Vent hole  
for c.f. NC /  
diaphragm sizes 10 - 50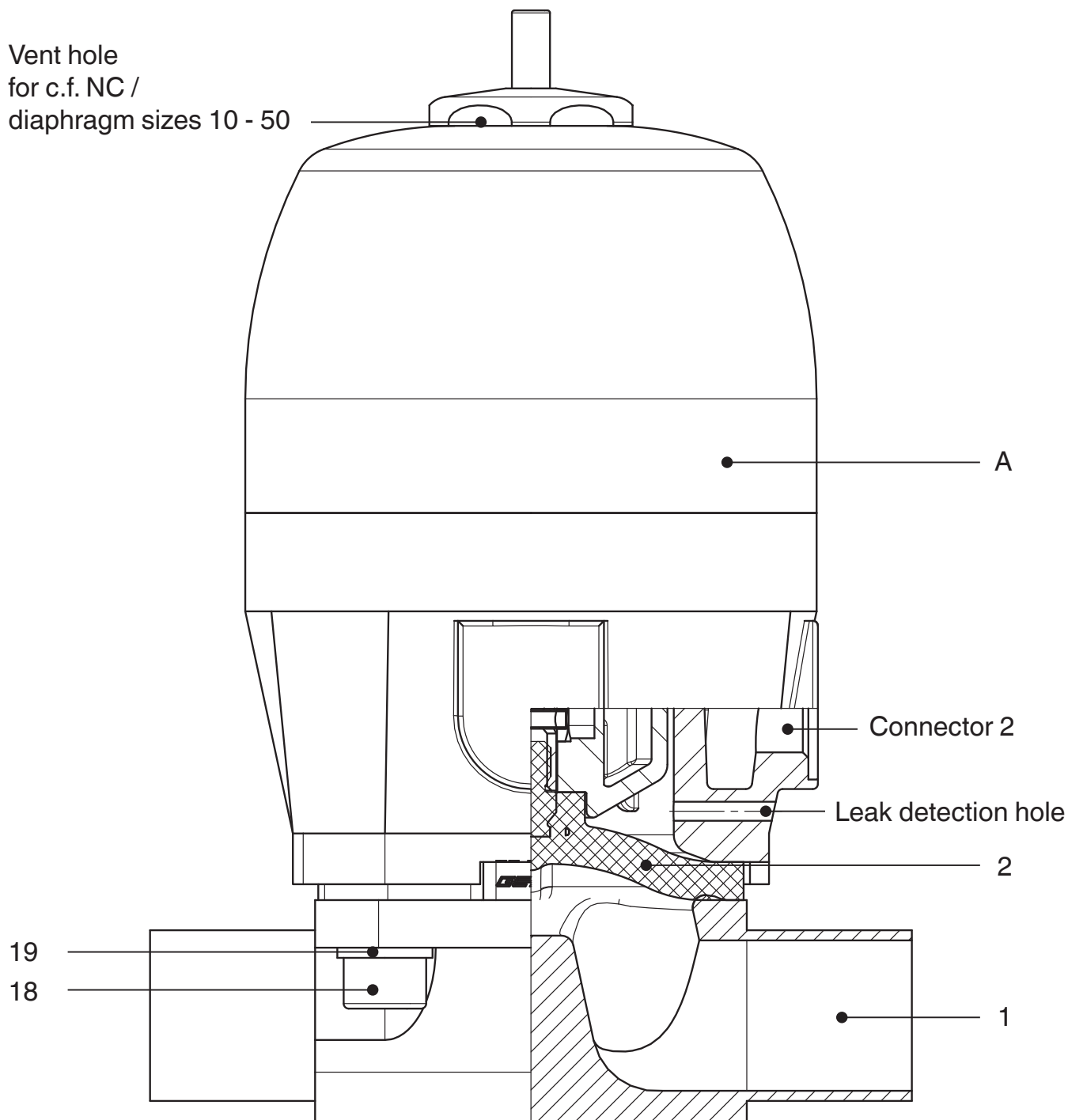

Item	Name	Order designation
Diaphragm sizes 10 - 50 / Actuator version: D		
1	Valve body	K600...
2	Diaphragm	600...M
18	Bolt	} 650...S30...
19	Washer	
A	Actuator	9650...

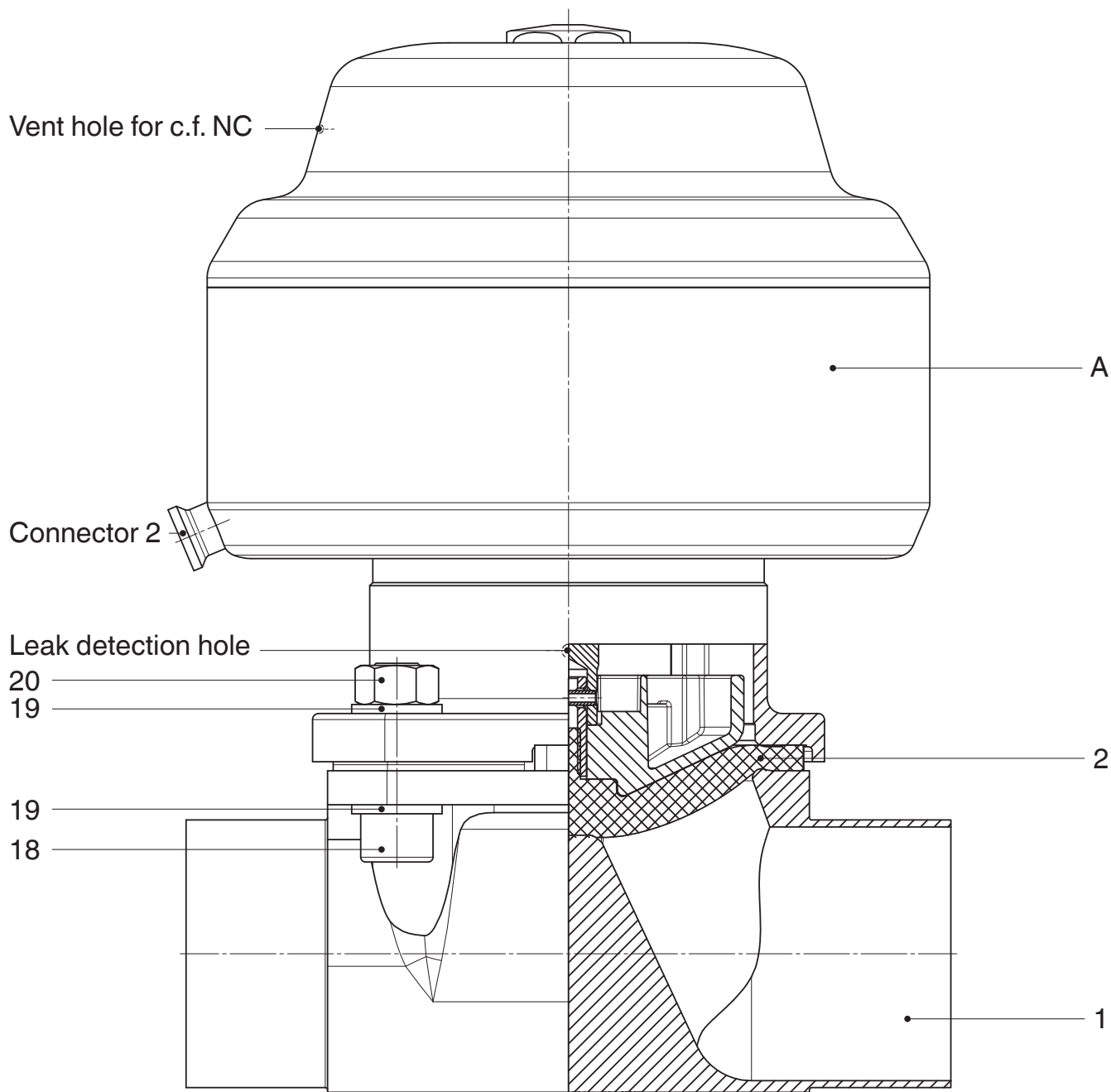Item	Name	Order designation
Diaphragm size 80		
1	Valve body	K600...
2	Diaphragm	600...M
18	Bolt	} 650...S30...
19	Washer	
20	Nut	
A	Actuator	9650...

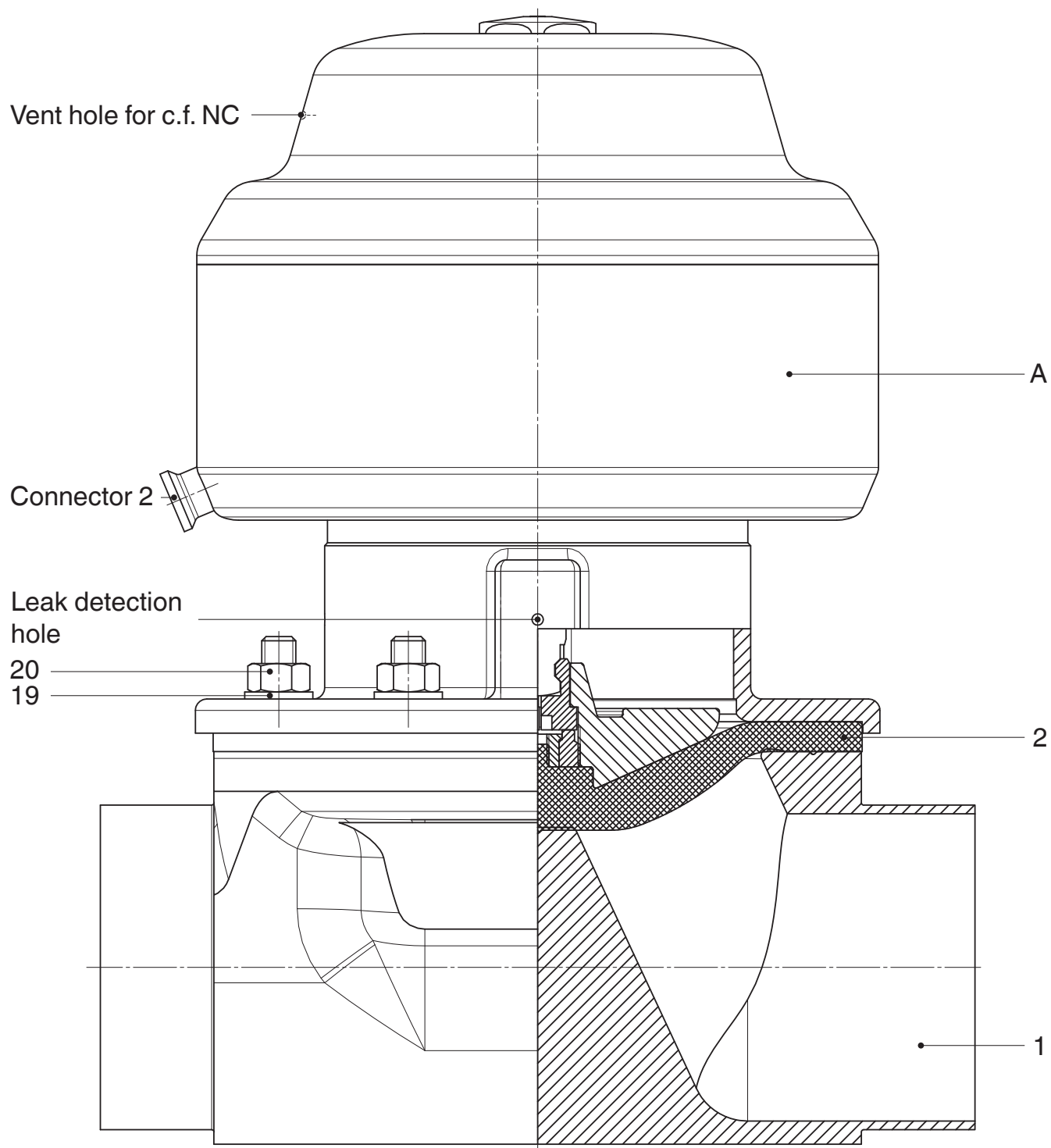Item	Name	Order designation
Diaphragm size 100		
1	Valve body	K600...
2	Diaphragm	600...M
19	Washer	} 650...S30...
20	Nut	
A	Actuator	9650...

# Declaration of Incorporation

**according to the EC Machinery Directive 2006/42/EC, Annex II, 1.B  
for partly completed machinery**

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**Description and identification of the partly completed machinery:**

Make: GEMÜ diaphragm valve, pneumatically operated  
Serial number: from 29.12.2009  
Project number: MV-Pneum-2009-12  
Commercial name: Type 650

**We hereby declare that the following essential requirements of the Machinery Directive 2006/42/EC have been fulfilled:**

1.1.3., 1.1.5., 1.2.1., 1.3., 1.3.2., 1.3.3., 1.3.4., 1.3.7., 1.3.9., 1.5.3., 1.5.5., 1.5.6., 1.5.7., 1.5.8., 1.5.9., 1.6.5.

**We also declare that the specific technical documentation has been compiled in accordance with part B of Annex VII.**

**We expressly declare that the partly completed machinery complies with the relevant provisions of the following EC directives:**


2006/42/EC:2006-05-17: (Machinery Directive) Directive 2006/42/EC of the European Parliament and of the Council of 17 May 2006 on machinery, and amending Directive 95/16/EC (recast) (1)

The manufacturer or his authorised representative undertake to transmit, in response to a reasoned request by the national authorities, relevant information on the partly completed machinery. This transmission takes place:

electronically

This does not affect the industrial property rights!

**Important note! The partly completed machinery may be put into service only if it was determined, where appropriate, that the machinery into which the partly completed machinery is to be installed meets the provisions of this Directive.**

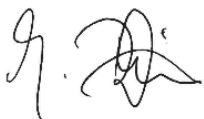Joachim Brien  
Head of Technical Department

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declare that the equipment listed below complies with the safety requirements of the Pressure Equipment Directive 2014/68/EU.

### Description of the equipment - product type

**Diaphragm valve**  
**GEMÜ 650**


Notified body:	TÜV Rheinland Industrie Service GmbH
Number:	0035
Certificate no.:	01 202 926/Q-02 0036
Technical standards used:	AD 2000

Conformity assessment procedure:  
**Module H1**

### Note for equipment with a nominal size $\leq$ DN 25:

The products are developed and produced according to GEMÜ process instructions and quality standards which comply with the requirements of ISO 9001 and ISO 14001.

According to Article 4, Paragraph 3 of the Pressure Equipment Directive 2014/68/EU, these products must not be identified by a CE-label.

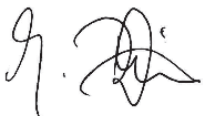Joachim Brien  
Head of Technical Department

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Hereby we, **GEMÜ Gebr. Müller Apparatebau GmbH & Co. KG**