

## GEMÜ D40

### Pneumatically operated diaphragm valve



#### Features

- Low-maintenance due to the firmly chambered diaphragm
- Fast, safe and simple diaphragm and actuator mounting
- Very high Kv values due to flow-optimised valve body
- Valve can be configured for the respective process parameters
- Identification of the angle of rotation (hash mark), optical position indicator and transparent cap as standard
- Simple modular expansion possible with future-orientated automation components
- Suitable for use in hygienic and aseptic applications (CIP/SIP capable and autoclavable)

#### Description

The pneumatically operated diaphragm valve GEMÜ D40 is designed for use in sterile applications. The diaphragm hermetically separates the actuator from the working medium. All actuator parts (except the seals and design elements) are made from stainless steel. The "Normally closed", "Normally open" and "Double acting" control functions are available.

#### Technical specifications

- **Media temperature :** -10 to 100 °C
- **Sterilization temperature:** max. 150 °C
- **Ambient temperature:** -10 to 80 °C
- **Operating pressure :** 0 to 16 bar
- **Nominal sizes:** DN 10 to 40
- **Body configurations:** Straight through body | T body
- **Connection types:** Clamp | Spigot
- **Connection standards :** ASME | DIN | EN | ISO
- **Body materials:** (316L), forging material | 1.4435 (316L), block material | 1.4435 (BN2), block material | 1.4435 (BN2), forged material
- **Diaphragm materials:** PTFE/EPDM
- **Conformities:** ATEX | FDA | Regulation (EC) No. 1935/2004 | USP

Technical data depends on the respective configuration




webcode: GM-D40



## Product description

### Construction

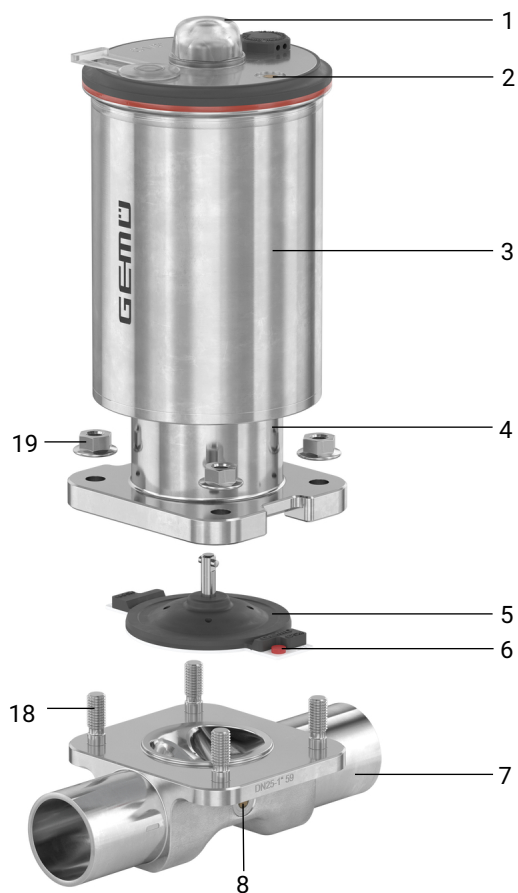

Position	Name	Materials
1	Transparent cap	PC
2	CONEXO actuator RFID chip (see Conexo information)	
3	Actuator	1.4301 / 1.4308
4	Distance piece with leak detection hole	1.4308 / 1.4408
5	Diaphragm	PTFE/EPDM two-piece
6	CONEXO diaphragm RFID chip (see Conexo information)	
7	Valve body	1.4435 (F316L), forged body 1.4435 (316L), block material 1.4435 (BN2), forged body, $\Delta$ Fe < 0.5% 1.4435 (BN2), block material, $\Delta$ Fe < 05 %
8	CONEXO body RFID chip (see Conexo information)	
18, 19	Screw connection kit (stud bolts, flange nuts)	

## GEMÜ CONEXO

The interaction of valve components that are equipped with RFID chips and an associated IT infrastructure actively increase process reliability.



The diaphragm is read out using an RFID reader, the CONEXO pen. The CONEXO app, which can be installed on mobile devices, not only facilitates and improves the “installation qualification” process, but also makes the maintenance process much more transparent and easier to document. The app actively guides the maintenance technician through the maintenance schedule and directly provides them with all the information assigned to the valve, such as test reports, testing documentation and maintenance histories. The CONEXO portal acts as a central element, helping to collect, manage and process all data.

**You can find further information on GEMÜ CONEXO at:**

[www.gemu-group.com/conexo](http://www.gemu-group.com/conexo)

## Range overviews

### Range overview of surface finishes

#### Internal surface finishes for block material body <sup>1)</sup>

Process contact surfaces	Mechanically machined <sup>2)</sup>		Electropolished	
	Hygiene class DIN 11866	Code	Hygiene class DIN 11866	Code
Ra ≤ 0.40 µm	H4	1536	HE4	1537

Process contact surfaces in accordance with ASME BPE <sup>3)</sup>	Mechanically machined <sup>2)</sup>		Electropolished	
	ASME BPE surface designation	Code	ASME BPE surface designation	Code
Ra max. = 0.51 µm (20 µinch)	SF1	SF1	-	-
Ra max. = 0.38 µm (15 µinch)	-	-	SF4	SF4

Ra acc. to DIN EN ISO 4288 and ASME B46.1

- 1) Surface finishes of customized valve bodies may be limited in special cases.
- 2) Or any other finishing method that meets the Ra value (acc. to ASME BPE).
- 3) When using these surfaces, the bodies are marked according to the specifications of ASME BPE.  
The surfaces are only available for valve bodies which are made of materials (e.g. GEMÜ material codes 40, 41) and use connections (e.g. GEMÜ connection codes 59, 80, 88) according to ASME BPE.

### Range overview of valve bodies

#### Range overviews of spigots 17, 59, 60

MG	DN	Connection type code <sup>1)</sup>		
		17	59	60
		Material code <sup>2)</sup>		
		40, 41, 42, 43		
<b>B</b>	<b>10</b>	X	-	X
	<b>15</b>	X	X	X
	<b>20</b>	X	X	-
	<b>25</b>	-	X	-
<b>C</b>	<b>20</b>	X	-	X
	<b>25</b>	X	X	X
	<b>32</b>	X	-	X
	<b>40</b>	X	X	-

MG = diaphragm size, X = standard

- 1) **Connection type, spigot 1**  
Code 17: Spigot EN 10357 series A / DIN 11866 series A formerly DIN 11850 series 2  
Code 59: Spigot ASME BPE/DIN 11866 series C  
Code 60: Spigot ISO 1127/EN 10357 series C/DIN 11866 series B
- 2) **Valve body material**  
Code 40: 1.4435 (F316L), forged body  
Code 41: 1.4435 (316L), block material  
Code 42: 1.4435 (BN2), forged body, Δ Fe < 0.5%  
Code 43: 1.4435 (BN2), block material, Δ Fe < 0.5%

**Range overviews of clamps 80, 82, 88, 8A, 8P, 8T**

MG	DN	Connection type code <sup>1)</sup>			
		80, 8P	82	88, 8T	8A
		Material code <sup>2)</sup>			
40, 41, 42, 43					
<b>B</b>	<b>10</b>	-	X	-	X
	<b>15</b>	X	X	X	X
	<b>20</b>	X	-	X	X
	<b>25</b>	X	-	X	-
<b>C</b>	<b>20</b>	-	-	-	-
	<b>25</b>	X	X	X	X
	<b>32</b>	-	X	-	X
	<b>40</b>	X	-	X	-

MG = diaphragm size, X = standard

1) **Connection type**

Code 80: Clamp ASME BPE, face-to-face dimension FTF ASME BPE, length only for body configurations D & F

Code 82: Clamp DIN 32676 Series B, Overall length FTF EN 558 Series 7, Overall length only for housing type D

Code 88: Clamp ASME BPE, for pipe ASME BPE, face-to-face dimension FTF EN 558 series 7, length only for body configurations D & F

Code 8A: Clamp DIN 32676 Series A, Overall length FTF according to EN 558 Series 7, Overall length only for housing type D

Code 8P: Clamp DIN 32676 series C, face-to-face dimension FTF ASME BPE, length only for body configurations D & F

Code 8T: Clamp DIN 32676 series C, face-to-face dimension FTF EN 558 series 7, length only for body configurations D & F

2) **Valve body material**

Code 40: 1.4435 (F316L), forged body

Code 41: 1.4435 (316L), block material

Code 42: 1.4435 (BN2), forged body,  $\Delta$  Fe < 0.5%

Code 43: 1.4435 (BN2), block material,  $\Delta$  Fe < 0.5%

## Order data

The order data provide an overview of standard configurations.

Please check the availability before ordering. Other configurations available on request.

## Order codes

1 Type	Code
Diaphragm valve, pneumatically operated, stainless steel piston actuator	D40

2 DN, connection 1	Code
DN 10	10
DN 15	15
DN 20	20
DN 25	25
DN 32	32
DN 40	40

3 Body configuration	Code
2/2-way body	D

4 Valve body connection type, connection 1	Code
Spigot EN 10357 series A/DIN 11866 series A	17
Spigot ASME BPE/DIN EN 10357 series C (from 2022 edition)/DIN 11866 series C	59
Spigot ISO 1127/DIN EN 10357 series C (2014 edition)/DIN 11866 series B	60
Clamp ASME BPE, for pipe ASME BPE, face-to-face dimension FTF ASME BPE, length only for body configuration D	80
Clamp DIN 32676, for pipe DIN 11866 series B, face-to-face dimension FTF EN 558 series 7, length only for body configuration D	82
Clamp ASME BPE, for pipe ASME BPE, face-to-face dimension FTF EN 558 series 7, length only for body configuration D	88
Clamp DIN 32676, for pipe DIN 11866 series A, face-to-face dimension FTF acc. to EN 558 series 7, length only for body configuration D	8A
Clamp DIN 32676, for pipe DIN 11866 series C, face-to-face dimension FTF ASME BPE, length only for body configuration D	8P
Clamp DIN 32676, for pipe DIN 11866 series C, face-to-face dimension FTF EN 558 series 7, length only for body configuration D	8T

5 Valve body material	Code
1.4435 (F316L), forged body	40
1.4435 (316L), block material	41
1.4435 (BN2), forged body, $\Delta$ Fe < 0.5%	42
1.4435 (BN2), block material, $\Delta$ Fe < 0.5%	43
Material code 41 and code 43 on request	

6 Diaphragm material	Code
PTFE/EPDM, two-piece	T1

7 Control function	Code
Normally closed (NC)	1
Normally open (NO)	2
Double acting (DA)	3

8 Actuator spring set	Code
Standard spring set	1

9 Diaphragm size	Code
Diaphragm size B	B
Diaphragm size C	C

10 Actuator size	Code
Actuator size 2	2
Actuator size 3	3
Actuator size 4	4

11 Valve body surface	Code
Ra ≤ 0.4 µm for media wetted surfaces, in accordance with DIN 11866 H4, mechanically polished internal mechanically machined internal	1536
Ra ≤ 0.4 µm for media wetted surfaces, in accordance with DIN 11866 HE4, electropolished internal/external	1537
Ra max. 0.51 µm (20 µin.) for media wetted surfaces, in accordance with ASME BPE SF1, mechanically machined internal	SF1
Ra max. 0.38 µm (15 µin.) for media wetted surfaces, in accordance with ASME BPE SF4, electropolished internal/external	SF4

12 CONEXO	Code
Integrated RFID chip for electronic identification and traceability	C
Without	

13 Special function	Code
ATEX marking	X

**Order example**

Ordering option	Code	Description
1 Type	D40	Diaphragm valve, pneumatically operated, stainless steel piston actuator
2 DN, connection 1	15	DN 15
3 Body configuration	D	2/2-way body
4 Valve body connection type, connection 1	17	Spigot EN 10357 series A/DIN 11866 series A
5 Valve body material	40	1.4435 (F316L), forged body
6 Diaphragm material	T1	PTFE/EPDM, two-piece
7 Control function	1	Normally closed (NC)
8 Actuator spring set	1	Standard spring set
9 Diaphragm size	B	Diaphragm size B
10 Actuator size	2	Actuator size 2
11 Valve body surface	1537	Ra ≤ 0.4 µm for media wetted surfaces, in accordance with DIN 11866 HE4, electropolished internal/external
12 CONEXO		Without
13 Special function	X	ATEX marking

## Technical data

### Medium

**Working medium:** Corrosive, inert, gaseous and liquid media which have no negative impact on the physical and chemical properties of the body and diaphragm material.

**Control medium:** Inert gases


### Temperature

Media temperature:	Diaphragm material	Standard
	PTFE/EPDM (code T1)	-10 to 100 °C

**Ambient temperature:** -10 – 80 °C

**Control medium temperature:** 0 – 60 °C

**Sterilization temperature:** PTFE/EPDM (code T1): Max. 150 °C, maximum 180 min per cycle  
 The sterilization temperature is valid for steam (saturated steam) or superheated water (max. 6 bar). PTFE diaphragms can also be used as moisture barriers; however, this will reduce their service life. It is not possible to retighten the connecting screws between the actuator and the valve body. This also applies to PTFE diaphragms exposed to high temperature fluctuations. The maintenance cycles must be adapted accordingly. The following valve arrangement has proven itself for interfaces between steam and process pipes: A globe valve for shutting off steam pipes and a diaphragm valve as an interface to the process pipes.

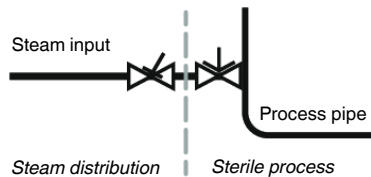

**Storage temperature:** -10 – 40 °C

### Pressure

**Operating pressure:** 0 – 16 bar

Information on operating pressures applied on both sides and for high purity media on request. Diaphragms with code T1 in valves in the 0–16 bar design have a reduced service life compared to the 0–10 bar design.

MG	DN	Actuator size	Diaphragm material	
			PTFE	
			Control function 1	Control function 2+3
<b>B</b>	<b>10 - 25</b>	<b>2</b>	0–10.0	0–10.0*
		<b>3</b>	0–16.0	-
<b>C</b>	<b>20 - 40</b>	<b>3</b>	0–5.0	0–5.0
		<b>4</b>	0–10.0	-

MG = diaphragm size

All pressures are gauge pressures. Operating pressure values were determined with static operating pressure applied on one side of a closed valve. Sealing at the valve seat and atmospheric sealing is ensured for the given values.

\* As a guide for operating the system with low wear on the diaphragm, refer to the control pressure – operating pressure diagrams below.

**Vacuum:** Can be used up to a vacuum of 70 mbar (absolute)

**Pressure rating:** PN 16

**Leakage rate:** Leakage rate A to P11/P12 EN 12266-1

**Filling volume:**

Actuator size	Diaphragm size	Control function		Valve stroke
		NC	NO	
2	B	0.06 dm <sup>3</sup>	0.05 dm <sup>3</sup>	7.5 mm
3	C	0.09 dm <sup>3</sup>	0.08 dm <sup>3</sup>	11.8 mm
4	C	0.181 dm <sup>3</sup>	0.193 dm <sup>3</sup>	11.8 mm

NC = filling volume when open, NO = filling volume when closed


**Control pressure:**

**Control function 1**

4.0 Up to 8.0 bar

**Control function 2 and 3: Control pressure/operating pressure diagram**


GEMÜ D40 design, diaphragm size B, actuator size 2

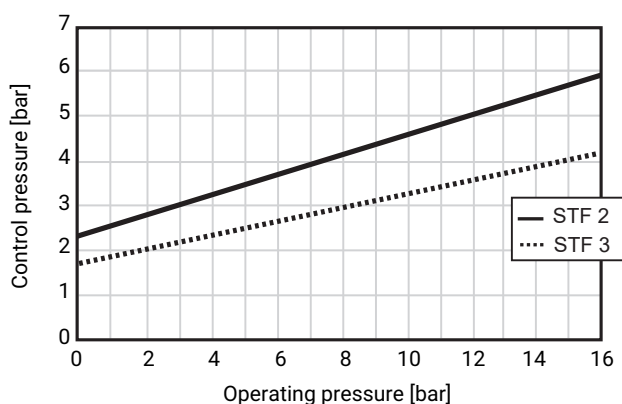

Stf. = control function

**Control function 2 and 3: Control pressure/operating pressure diagram**

GEMÜ D40 design, diaphragm size C, actuator size 4

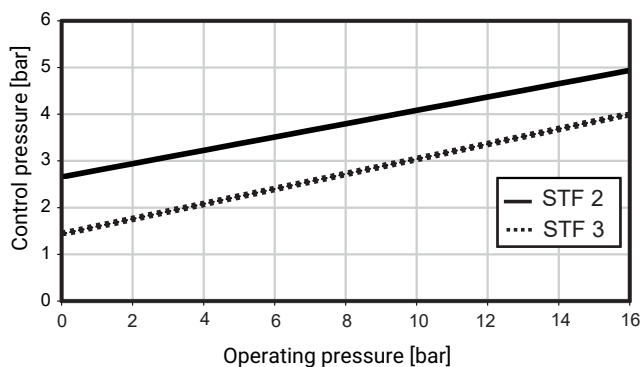

Stf. = control function

The control pressure depending on the prevailing operating pressure, as shown in the diagram, is intended as a guide for operating the system with low wear on the diaphragm.

**Kv values:**

Diaphragm size	DN	Connection type (code)		
		17	59	60
<b>B</b>	<b>10</b>	4.4	-	8.0
	<b>15</b>	9.5	3.9	9.5
	<b>20</b>	9.5	9.5	-
	<b>25</b>	-	9.5	-
<b>C</b>	<b>20</b>	16.5	-	23.2
	<b>25</b>	25.3	22.0	26.4
	<b>32</b>	26.4	-	24.3
	<b>40</b>	24.6	25.2	-

MG = diaphragm size, Kv values in m³/h


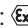
Kv values determined in accordance with EN 60534, inlet pressure 5 bar, Δp 1 bar, stainless steel valve body and two-piece PTFE/EPDM diaphragm. The Kv values for other product configurations (e.g. other diaphragm or body materials) may differ. In general, all diaphragms are subject to the influences of pressure, temperature and process parameters, as a result of which the Kv values deviate beyond the tolerance limit of the standard. The Kv value curve (Kv value dependent on valve stroke) can vary depending on the diaphragm material and term of use.

**Product conformities**

**Machinery Directive:** 2006/42/EC

**Pressure Equipment Directive:** 2014/68/EU

**Explosion protection:** ATEX (2014/34/EU), order code Special version X

**ATEX marking:** Gas:  II 2 G Ex h IIC T6 ... T3 Gb X  
Dust :  II -/2 D Ex h -/IIIC T150 °C -/Db X

**Food:** FDA  
USP Class VI  
Regulation (EC) No. 1935/2004 (for material code 40 only)  
Regulation (EC) No. 10/2011

**Mechanical data**

**Weight:**

**Body**

MG	DN	Spigot		
		Connection type code		
		17	59	60
<b>B</b>	<b>10</b>	0.33	-	0.32
	<b>15</b>	0.31	0.34	0.3
	<b>20</b>	0.28	0.31	-
	<b>25</b>	-	0.27	-
<b>C</b>	<b>20</b>	1.00	-	0.97
	<b>25</b>	0.94	0.98	0.89
	<b>32</b>	0.86	-	0.80
	<b>40</b>	0.78	0.81	-

Weights in kg  
MG = diaphragm size

**Weight:**

**Actuator (control function 1, NC)**


MG	Actuator size	Weight
<b>B</b>	2	1.7
<b>B</b>	3	1.2
<b>C</b>	3	1.4
<b>C</b>	4	3.8

Weights in kg  
MG = diaphragm size

**Angle of rotation:**


MG	DN	Connection type (code)		
		17	59	60
		$\alpha$		
<b>B</b>	<b>10</b>	36.3°	-	23.4°
	<b>15</b>	18.0°	38.1°	12.3°
	<b>20</b>	7.9°	18.3°	-
	<b>25</b>	-	3.0°	-
<b>C</b>	<b>20</b>	35.4°	-	28.4°
	<b>25</b>	24.5°	31.2°	18.1°
	<b>32</b>	14.9°	-	5.9°
	<b>40</b>	7.9°	10.8°	-

MG = diaphragm size

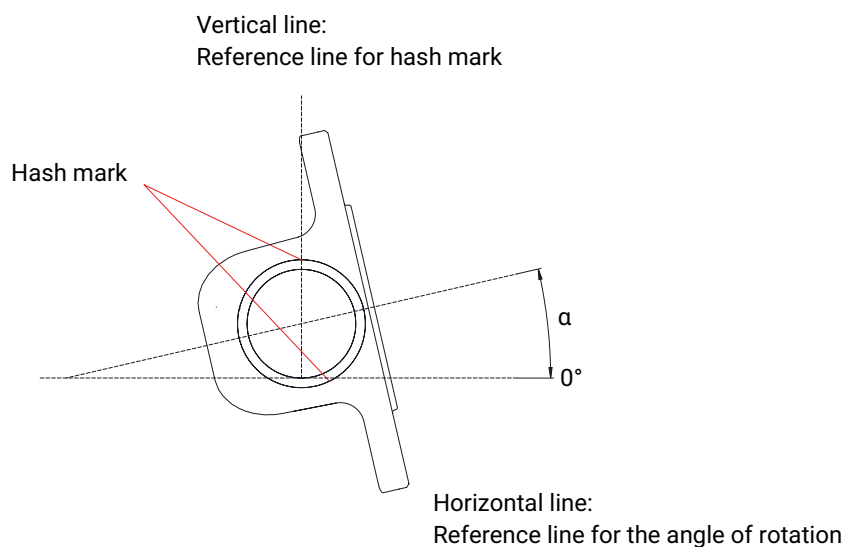

## Dimensions

### Actuator dimensions

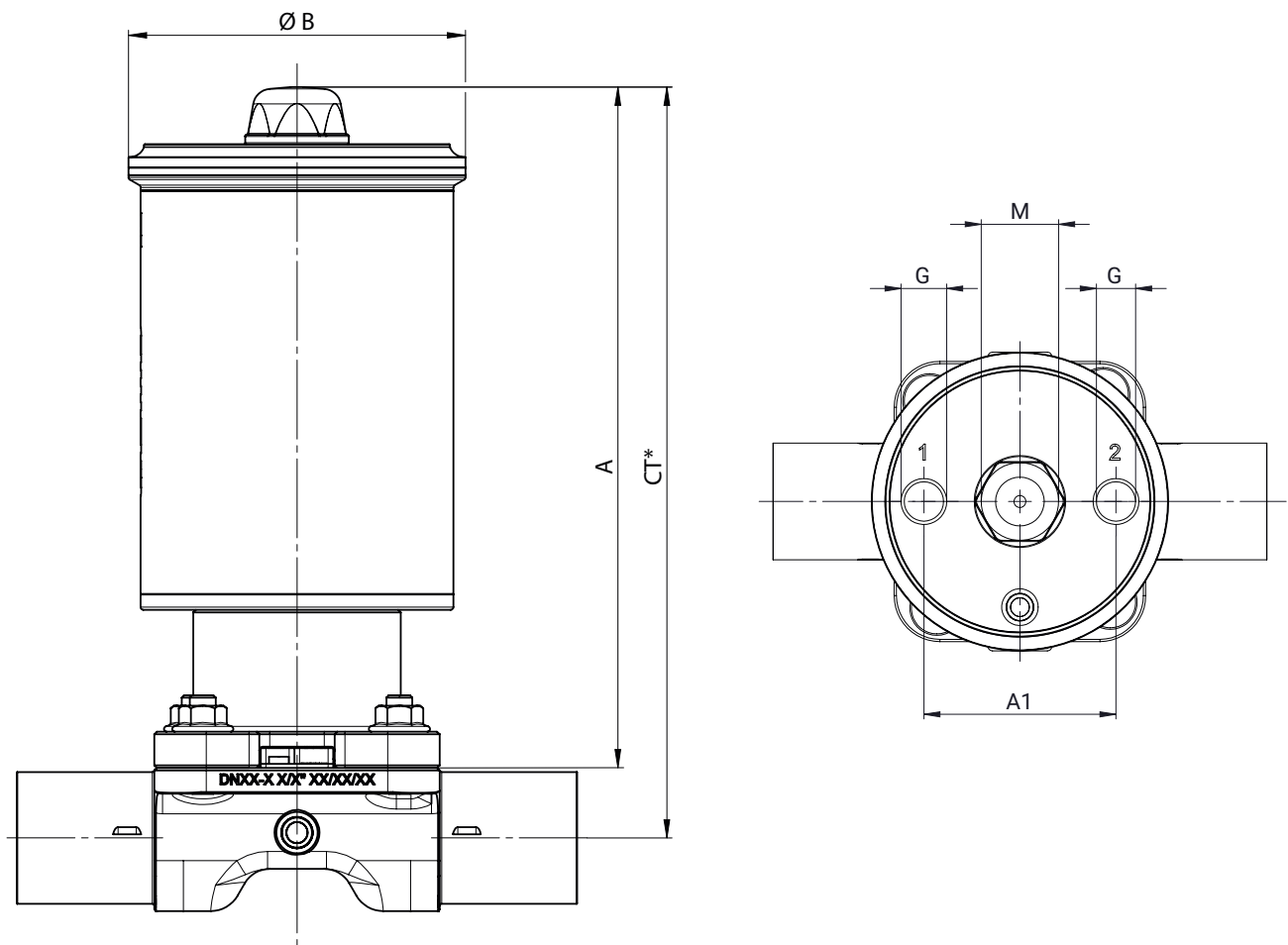

Actuator size	MG	A	$\varnothing B$	A1	G	M
2	B	131.3	65.0	42.0	G 1/8	M16x1
3	B	136.8	70.0	42.0	G 1/8	M16x1
3	C	154.1	70.0	42.0	G 1/8	M16x1
4	C	186.5	88.9	61.0	G 1/4	M26x1.5

Dimensions in mm


MG = diaphragm size

\* CT = A + H1 (see body dimensions)

## Body dimensions

### Spigot DIN/EN/ISO/ASME (code 17, 59, 60), forged material (code 40, 42)

Connection type spigot DIN/EN/ISO/ASME (codes 17, 59, 60)<sup>1)</sup>, forged material (code 40)<sup>2)</sup>

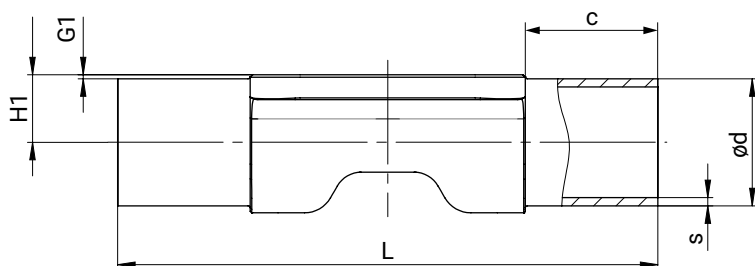

MG	DN	Pipe standard			EN 10357 series A/DIN 11866 series A	ASME BPE/DIN 11866 series C	ISO 1127/EN 10357 series C/DIN 11866 series B						
					Connection code								
					17	59	60						
		L	c (min)	G1	H1	Ød	s	H1	Ød	s	H1	Ød	s
<b>B</b>	<b>10</b>	108.0	25.0	0.8	7.3	13.0	1.5	5.57	9.53	0.89	9.4	17.2	1.6
	<b>15</b>	108.0	25.0	0.8	10.3	19.0	1.5	7.15	12.70	1.65	11.45	21.3	1.6
	<b>20</b>	108.0	25.0	0.8	12.3	23.0	1.5	10.33	19.05	1.65	-	-	-
	<b>25</b>	108.0	25.0	0.8	-	-	-	13.50	25.40	1.65	-	-	-
<b>C</b>	<b>20</b>	127.0	25.0	0.8	12.3	23.0	1.5	-	-	-	14.25	26.9	1.6
	<b>25</b>	127.0	25.0	0.8	15.3	29.0	1.5	13.5	25.4	1.65	17.65	33.7	2.0
	<b>32</b>	127.0	25.0	0.8	18.3	35.0	1.5	-	-	-	22.0	42.4	2.0
	<b>40</b>	127.0	25.0	0.8	21.3	41.0	1.5	19.85	38.1	1.65	-	-	-

Dimensions in mm

MG = diaphragm size

#### 1) Connection type

Code 17: Spigot EN 10357 series A/DIN 11866 series A, formerly DIN 11850 series 2

Code 59: Spigot ASME BPE/DIN EN 10357 series C (from 2022 edition)/DIN 11866 series C

Code 60: Spigot ISO 1127/DIN EN 10357 series C (2014 edition)/DIN 11866 series B


#### 2) Valve body material

Code 40: 1.4435 (F316L), forged body

Code 42: 1.4435 (BN2), forged body,  $\Delta Fe < 0.5\%$

**Clamp DIN/EN/ISO/ASME (codes 80, 82, 88, 8A, 8P, 8T), forged material (code 40, 42)**

Connection type clamp DIN/ASME (codes 80, 88, 8P, 8T)<sup>1)</sup>, forged material (code 40, 42)<sup>2)</sup>

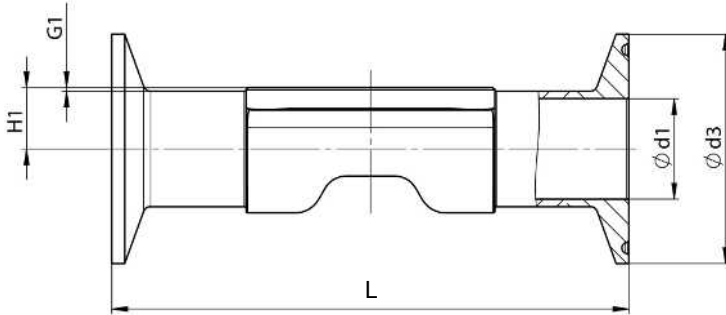

MG	DN	NPS	ød1		ød3		H1	G1	L	
			Connection type		Connection type				Connection type	
			80, 8P	88, 8T	80, 8P	88, 8T			80, 8P	88, 8T
<b>B</b>	<b>10</b>	3/8"	-	-	-	-	-	0.8	-	-
	<b>15</b>	1/2"	9.40	9.40	25.00	25.00	7.15	0.8	88.90	108.00
	<b>20</b>	3/4"	15.75	15.75	25.00	25.00	10.33	0.8	101.60	117.00
	<b>25</b>	1"	22.10	22.10	50.50	50.50	13.50	0.8	114.30	127.00
<b>C</b>	<b>20</b>	3/4"	-	-	-	-	-	0.8	-	-
	<b>25</b>	1"	22.10	22.10	50.50	50.50	13.50	0.8	114.30	127.00
	<b>32</b>	1 1/4"	-	-	-	-	-	0.8	-	-
	<b>40</b>	1 1/2"	34.80	34.80	50.50	50.50	19.85	0.8	139.70	159.00

Dimensions in mm

MG = diaphragm size

**1) Connection type**

Code 80: Clamp ASME BPE, face-to-face dimension FTF ASME BPE, length only for body configuration D

Code 88: Clamp ASME BPE, for pipe ASME BPE, face-to-face dimension FTF EN 558 series 7, length only for body configuration D

Code 8P: Clamp DIN 32676 series C, face-to-face dimension FTF ASME BPE, length only for body configuration D

Code 8T: Clamp DIN 32676 series C, face-to-face dimension FTF EN 558 series 7, length only for body configuration D

**2) Valve body material**

Code 40: 1.4435 (F316L), forged body

Code 42: 1.4435 (BN2), forged body, Δ Fe < 0.5%

**Connection type clamp DIN/ISO (code 82)<sup>1)</sup>, forged material (code 40, 42)<sup>2)</sup>**

MG	DN	NPS	ød1	ød3	H1	G1	L
			Connection type	Connection type			Connection type
			82	82			82
<b>B</b>	<b>10</b>	3/8"	14.00	25.00	9.4	0.8	108.00
	<b>15</b>	1/2"	18.10	50.50	11.45	0.8	108.00
	<b>20</b>	3/4"	-	-	-	0.8	-
	<b>25</b>	1"	-	-	-	0.8	-
<b>C</b>	<b>20</b>	3/4"	23.00	50.50	14.25	0.8	117.00
	<b>25</b>	1"	29.70	50.50	17.65	0.8	127.00
	<b>32</b>	1¼"	38.40	64.00	22	0.8	146.00
	<b>40</b>	1½"	-	-	-	0.8	-

Dimensions in mm

MG = diaphragm size

**1) Connection type**

Code 82: Clamp DIN 32676 Series B, Overall length FTF EN 558 Series 7, Overall length only for housing type D

**2) Valve body material**

Code 40: 1.4435 (F316L), forged body

Code 42: 1.4435 (BN2), forged body, Δ Fe &lt; 0.5%

**Connection type clamp DIN/EN (code 8A)<sup>1)</sup>, forged material (code 40, 42)<sup>2)</sup>**

MG	DN	NPS	ød1	ød3	H1	G1	L
			Connection type	Connection type			Connection type
			8A	8A			8A
<b>B</b>	<b>10</b>	3/8"	10.00	34.00	7.30	0.8	108.00
	<b>15</b>	1/2"	16.00	34.00	10.30	0.8	108.00
	<b>20</b>	3/4"	20.00	34.00	12.30	0.8	117.00
	<b>25</b>	1"	-	-	-	0.8	-
<b>C</b>	<b>20</b>	3/4"	20.00	34.00	12.30	0.8	117.00
	<b>25</b>	1"	26.00	50.50	15.30	0.8	127.00
	<b>32</b>	1¼"	32.00	50.50	18.30	0.8	146.00
	<b>40</b>	1½"	38.00	50.50	21.30	0.8	159.00

Dimensions in mm

MG = diaphragm size

**1) Connection type**

Code 8A: Clamp DIN 32676 Series A, Overall length FTF according to EN 558 Series 7, Overall length only for housing type D

**2) Valve body material**

Code 40: 1.4435 (F316L), forged body

Code 42: 1.4435 (BN2), forged body, Δ Fe &lt; 0.5%

## Accessories



### GEMÜ 12A0

#### Intelligent electrical position indicator

Independent of the actuator size and control function, the GEMÜ 12A0 electrical position indicator, as an automation module, is compatible with all pneumatically operated process valves of the new valve generation and with quarter turn valves. Contactless position detection determines the valve position precisely, reliably and without being subject to wear. The current valve position is displayed via high visibility LEDs, and fed back via electrical signals. In addition to this, there is an integrated mechanical position indicator. Modern communication interfaces, an integrated sensor system and the GEMÜ app operating option are all features that characterize the innovative electrical position indicator.



### GEMÜ 44A0

#### Multi-functional valve actuation

Independent of the actuator size, the GEMÜ 44A0 multi-functional valve actuation, as an automation module, is compatible with all pneumatically operated process valves with single acting linear actuator of the new valve generation. Depending on the order variant and the set device functions, the connected process valves can be controlled conventionally open/closed (combi switchbox) or the valve position can be precisely controlled (positioner). Contactless position detection determines the valve position precisely, reliably and without being subject to wear. The current valve position is displayed via high visibility LEDs, and fed back via electrical signals. In addition to this, there is an integrated mechanical position indicator. Modern communication interfaces, an integrated sensor system and the GEMÜ app operating option are all features that characterize this innovative product.



GEMÜ Gebr. Müller Apparatebau GmbH & Co. KG  
Gert-Müller-Platz 1, 74635 Kupferzell, Germany  
Phone +49 (0) 7940 1230 · info@gemue.de  
www.gemu-group.com